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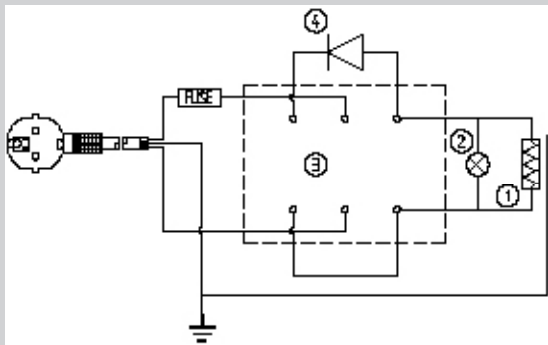
CAUTION

Read these instructions carefully before installing and starting up the equipment.

Do not touch the heating element. Burns can be caused due to its high temperature.

Do not use the equipment if it is not earthed.

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This product is guaranteed for one year. The guarantees does not cover damages caused by a misuse of the equipment or for reasons beyond J.P. SELECTA, s.a.u.

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ACCESSORIES

Stand bar, in Duraluminium, Ø 12mm. 700mm length.
 Code: 6000270

**MANTAS CALEFACTORAS PARA MATRACES****HEATING MANTLES FOR FLASKS****“Fibroman-C”**

3003141 3003142
 3003145 3031410
 3031420 3031430
 3031440 3031450
 3031460 3314100
 3314200*

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*The mantle part no 3314200 is supplied without the 2 position controller and the switch.

Código/ Code	Capacidad /Capacity ml	Potencia /Power W	Peso/Weight (Kg.)
3003141	100	125	0.8
3003142	250	120	0.9
3003145	500	250	1
3031410	1000	475	1.2
3031420	2000	700	1.8
3031430	3000	750	2
3031450	5000	1500	3
3314100	10000	2250	4.5
3314200	20000	2800	11

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FEATURES

- *Temperature up to 450°C on mantle surface.*
- *For spheric flasks from 100ml to 20.000ml.*
- *Woven mineral glass fiber, fully free of asbestos.*
- *Heating element evenly distributed through the body of the mantle.*
- *Two power selector.*

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Notice to customers:

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.

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INSTRUCTIONS

- *During the first starting up, you will observe that the heating mantle gives off some fumes and a characteristic smell; it is due to the dressing of the heating material. These circumstances are completely normal and will disappear in a period of time.*
- *It is recommended to use adjustable flasks with the mantle.*
- *Avoid spilling liquids on the heating material surface, since it means the immediate damage of the mantle.*
- *Make sure that the mains power corresponds to the one indicated on the description plate.*
- *For your safety, do not switch on the apparatus if it is not earthed.*
- *Turn on the main switch «POWER» (1) on selecting the desired power (50% or 100%).*