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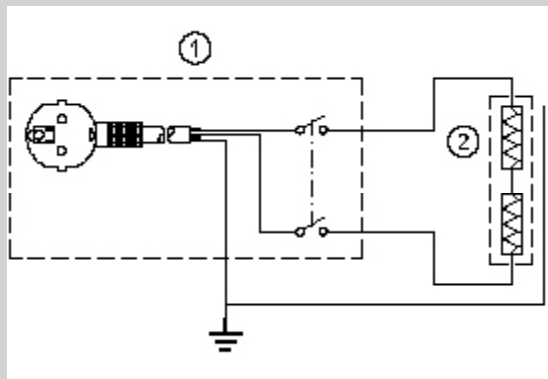
CAUTION

Read these instructions carefully before installing and starting up the equipment.

Do not touch the heating element. Burns can be caused due to its high temperature.

Do not use the equipment if it is not earthed.

COMPONENTES DE RECAMBIO / SPARE PARTS

DESCRIPCIÓN / DESCRIPTION:	COD.
Cable + selector	7070
MANTA / Heating mantle 250 ml.	41001
MANTA / Heating mantle 500ml.	41002
MANTA / Heating mantle 1000ml.	41003
MANTA / Heating mantle 2000ml.	41004

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This product is guaranteed for one year. The guarantees does not cover damages caused by a misuse of the equipment or for reasons beyond J.P. SELECTA, s.a.u.

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FEATURES

- Temperature up to 450°C on the mantle surface.
- For spheric flasks from 250 to 2.000ml.
- Woven mineral glass fiber, fully free of asbestos.
- Heating element evenly distributed throughout the body of the mantle.
- With attachment for ring or tripod stand.

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Notice to customers:

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.

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INSTRUCTIONS

- During the first starting up, you will observe that the heating mantle gives off some fumes and a characteristic smell; it is due to the dressing of the heating material. These circumstances are completely normal and will disappear in a period of time.
- It is recommended to use adjustable flasks with the mantle.
- Avoid spilling liquids on the heating material surface, since it means the immediate damage of the mantle.
- Make sure that the mains power corresponds to the one indicated on the description plate.
- For your safety, do not switch on the apparatus if it is not earthed.
- Turn on the main switch «POWER» (1).