DISPENSER

4000490

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the apparatus with liquids which can give off vapours capable of making explosive mixtures.*

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The standard equipment consist of the following components:

| Pieza | Piece | Código / Code | Cant. / Qty. |
|-------------------------|--------------------|---------------|--------------|
| Dispenser | Dispenser | 4000490 | 1 |
| Tapa | Cover | 30012 | 1 |
| Manual de instrucciones | Instruction manual | 80051 | 1 |
| Termómetro | Thermometer | 1000100 | 1 |

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Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

| Código / Code | Capacidad litros / Capacity litres | Medidas útiles / Tank dimensions cm | | Medidas exteriores / Overall dimensions cm | | Consumo / Consump. (W) | Estabilidad * / Stability * (K) | Peso / Weight (Kg) |
|---------------|------------------------------------|-------------------------------------|--------------|--|---------------|------------------------|---------------------------------|--------------------|
| | | Ø | Alto / Depth | Ø | Alto / Height | | | |
| 4000490 | 4 | 20 | 15 | 28 | 30 | 400 | ±1.5 | 5 |

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* Measured according to the measurement procedure of J.P. SELECTA, s.a. u.

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Paraffin wax DISPENSER are designed for adjustable temperatures from ambient +5°C up to 100°C (Stability ±1.5°C).

Epoxy-coated outer casing, inner tank and lid with PTFE-coated aluminium indelible matt black.

Tap dispenser with built-in heater.

Control panel:

- Main switch with indicator lamp.
- Heater operation indicator lamp.
- Temperature regulator thermostat.
- Tap temperature regulator.

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INSTALLATION

Place paraffin wax on a flat horizontal, level surface, leaving a space at the back and sides of the apparatus of about 10cm.

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Be sure that the voltage supply is the same as the one indicated on the characteristics plate.

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Do not use the autoclave if it is not earthed.

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If you change the plug bear in mind the following:

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Blue cable: Neutral.
Brown cable: Phase.
Yellow/green cable: Earth.

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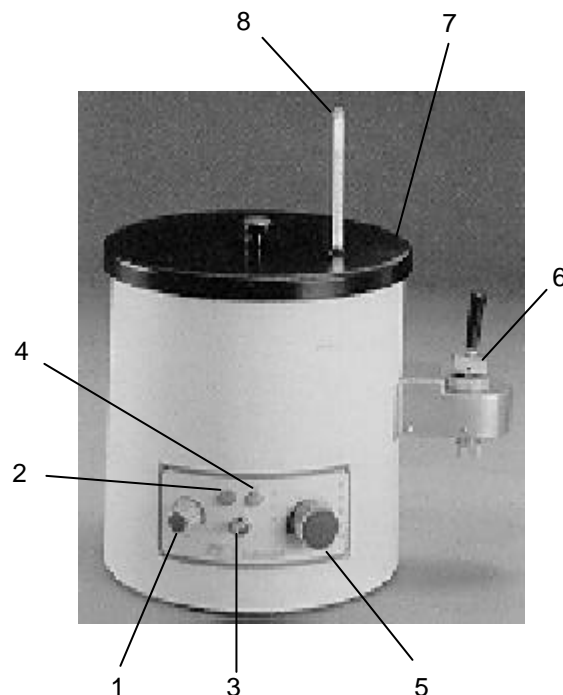
STARTING UP:

1. *Plug in the bath to the socket, with the main switch and the thermostat knob in "0" position.*
2. *Fill the bath with paraffin pills up to the desired level, bearing in mind that when is liquid, the paraffin reduces.*
3. *Turn the main switch (3) to the position I, at the same time the mains indicator lamp (green) (2) will light up.*
4. *Turn the temperature selector (5) to the desired temperature. At the same time the heater operation indicator lamp (yellow) (4) will light up.*
5. *After a certain time, the heater operation indicator lamp (4), will start operating intermittently. When these intermittences become uniform, the temperature is stabilized.*
6. *Although the thermostats controls the selected temperature, it will observe that the paraffin smelts slowly from bottom to up, so that up to it is smelted completely the temperature of the bath cannot be read with a thermometer. Once the reading of the complete liquid paraffin is done, if it is not the desired one, turn the thermostat knob left or right as required, and let it stabilize again.*
7. *The thermostat scale is intended as a guide, the real temperature level is the one that will show a thermometer introduced inside the bath.*
8. *The bath has to be always filled of paraffin. Replenish when necessary.*

REGULATION OF THE OUTLET TAP TEMPERATURE:

The regulation is done by turning the knob (1). Depending on the paraffin temperature, this knob will be moved. At first, it will be placed at maximum, and once the paraffin has liquated and it leaks out of the tap, the temperature can be reduced. Should the paraffin is solidified in the tap, the temperature must be increased again.

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MAINTENANCE

Before removing the cover disconnect the apparatus from the mains.

The manipulation of the internal electronic circuits of the apparatus by unauthorized personnel can cause irreparable damage. Take it to one of the J.P.SELECTA authorized technical services.

CLEANING:

For the cleaning of the different parts use the following products:

Cleaning of stainless steel: Alcohol

Cleaning of plastic: Alcohol with cotton duster.

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SPARE PARTS

To guarantee the safety of the equipment, the spare parts must be purchased from J.P.SELECTA,S.A.

| Descripción | Description | Código / Code |
|--------------------------|----------------------|---------------|
| Modelo | Model | 4000490 |
| Botón mando | Control knob | 6006 |
| Tapa botón mando | Cover knob | 6007 |
| Flecha botón mando | Arrow knob | 6008 |
| Alargo Schuko | Schuko cable | 7002 |
| Resistencia grifo | Tap heating element | 39077 |
| Resistencia baño | Bath heating element | 39082 |
| Termostato de regulación | Thermostat | 43006 |

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GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the apparatus by personnel not authorized by J.P. SELECTA, s.a.u. automatically cancels the guarantee.

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