**ULTRASONS-P**

**3000910**

**BAÑO LIMPIADOR DE PIPETAS  
PIPETTES ULTRASONIC CLEANER**

## INFORMACIÓN GENERAL

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## GENERAL INFORMATION

*Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*

*Read this handbook before use.*

*If you have any doubts, contact with J.P. Selecta.*

*If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the usage directive. The manufacturer is not responsible for the damage that can occur.*

*Do not use the bath with liquids which can give off vapours capable of making explosive mixtures.*

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## PACKING LIST

*The standard equipment consist of the following components:*

Descripción	Description	Código / Code	Cant. / Qty.
ULTRASONS-P	ULTRASONS-P	3000910	1
Manguera P.V.C.	P.V.C. tube	46080	2
Brida Mikalor	Brida Mikalor	44288	4
Canasta portapipeta	Basket with handle	28088	1
Cable de conexión	Connection cable	7117	1
Tapa	Cover	14100	1
Manual de instrucciones	Instruction manual	80088	1

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## ACCESSORIES

\*Detergent **MICRO**

*Specially for cleaning with **ULTRASONS** ultrasonic cleaners. We advise of this special detergent by its excellent cleaning power, it is biodegradable and free of phosphates. It is very useful for cleaning electronic components, pieces of glass, metal, plastic, rubber, jewels and materials with hard surfaces in general.*

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*Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.*

Código / Code	Capacidad litros / Capacity litres	Medidas útiles cm / Cylindrical vessel size cm		Medidas exteriores cm / Overall dimensions cm			Potencia W. / Generator power W	Reloj / Timer	Antiparásito / Antiparasite	Peso Kg. / Weight Kg
		Alto / Depth	Ancho / Length	Alto / Height	Ancho / Length	Fondo / Width				
3000910	10	64	16Ø	84	34	40	250	0-15'	Si / Yes	15

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## INSTALLATION

- *Place the appliance on a flat surface or on the ground, near a water intake and a sump. Leave a free space of 30cm around the equipment.*
- *Connect the water inlet in (5).*
- *Connect the outlet to the drain in (4) fixed with flanges.*
- *Ensure that the outlet does not form pockets or elbows that hinder the free circulation of water.*
- *Do not start the equipment without water.*

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*Do not use inflammable solvents or which tends to form explosive gases.*

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*Never leave the apparatus on without liquid in the tank as it could cause breakdowns.*

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*Be sure that the voltage supply is the same as the one indicated on the characteristics plate.*

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*Do not use the apparatus if it is not earthed.*

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*If you change the plug bear in mind the following:*

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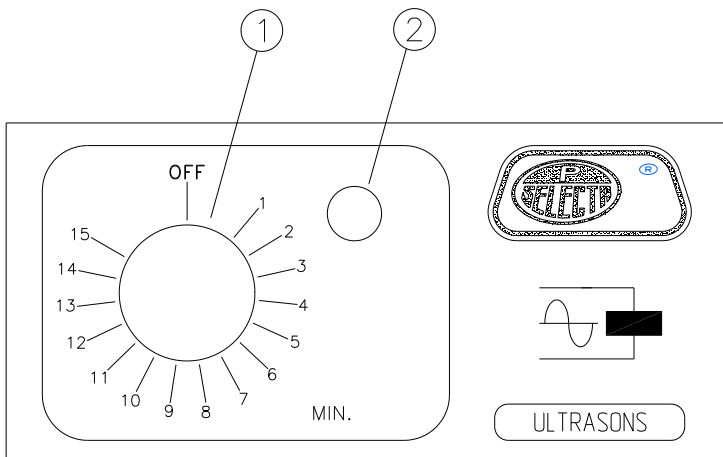
**Blue cable: Neutral  
Brown cable: Phase  
Yellow/green cable: Earth**

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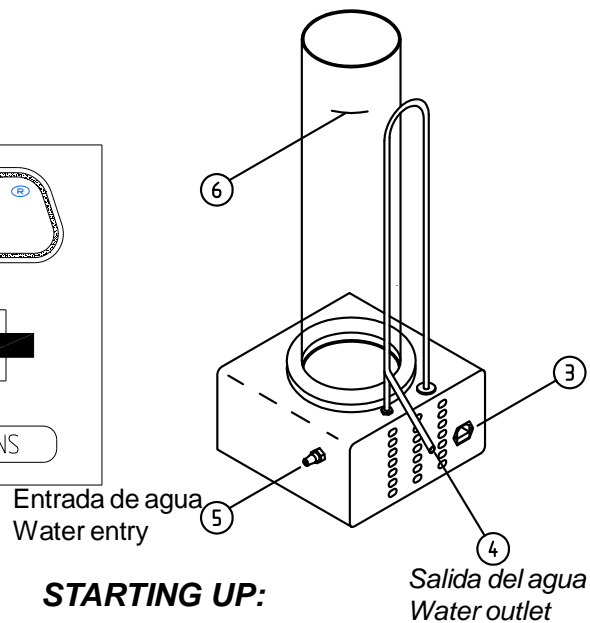
### STARTING UP:

1. Place the pipettes in vertical position inside the basket.
2. Place the basket with the pipettes inside the bath tank.
3. Turn on the water tap until the level (6) is reached. Additives may be added during the filling to facilitate cleaning.
4. Start up by using the timer knob (1) clockwise for temporasing mode.
5. At the desired time, open the water supply tap and adjust it so that the water level is maintained above siphon curve. This operation primes the siphon producing a circulation which causes the emptying of the cleaning liquid and constant renewal of the water (rinsed).
6. Once the pipettes are rinsed, turn off the apparatus by moving the timer knob (1) to "OFF" and turning off the water tap. Due to the siphon effect, the tank will be completely emptied and the apparatus will be ready for re-use.

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**COMMENTS:**

- \* *Do not overload the cleaner.*
- \* *Avoid placing heavy instruments directly on the bottom of tank (use the baskets provided for this purpose).*
- \* *Before cleaning anything -particularly valuable items- for the first time, it is worth experimenting with a sole piece before continuing with the others. Organic pearls are unsuitable for ultrasonic cleaners. Heat is generated during cleaning which could harm them.*
- \* *The chemical cleaning solutions employed must be compatible with work tools and furniture.*
- \* *Corrosive solvents must be avoided.*
- \* *usually 5-15 minutes is suitable time to clean.*
- \* *For greater conservation and durability of cleaner, it is better to keep it connected only during the cleaning of items.*
- \* Use demineralized water.

**MAINTENANCE**

**Before removing the casing, disconnect the apparatus from the mains.**

**The control panel must only be manipulated by authorized personnel.**

**GUARANTEE**

*This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use of causes beyond the control of J.P. SELECTA,S.A..*

*Any manipulation of the apparatus by unauthorized personnel by J.P. SELECTA,S.A. cancels the guarantee automatically.*

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To guarantee the safety of the equipment, the spare parts must be purchased from J.P.SELECTA,S.A.**