LABOR-LAM

3000438

**LÁMPARA DE LABORATORIO
LABORATORY LAMP**

INFORMACIÓN GENERAL

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.*

PACKING LIST

The standard equipment consists of the following components:

Piece	Code	Qty.
<i>Labor-Lam</i>	3000438	1
<i>Instruction manual</i>	80103	1

ACCESSORIES

STAND TO USE IN SLANTING POSITION, for examination of films, transparencies, analysis, etc...

For Labor-lam 3000438

Code 7000483

<i>Stainless steel hinge screw</i>	2 units
<i>Aluminium separator</i>	2 units
<i>Fixing wing nut</i>	2 units

TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz.

Code	Screen useful measures cm		Overall dimensions cm			Sreen max. temp.	Max. inclination	LED lamp	Weight Kg.
	Height	Length	Height	Length	Width				
3000438	8.5	30	32	13	5.5	35°C	45°C	12V 6000K	1.4

EQUIPMENT DESCRIPTION

LABOR-LAM lamps are recommended for film examinations, transparencies, graphs, analysis, etc.
Equipped with a clip stand for transparencies, plastic screen with rounded edges to avoid scratches.

INSTALLATION AND STARTING UP

* Installation in vertical position:

- Fix the lamp onto the holes at the back of the casing.
- Connect the plug into the mains.
- Press the switch to start the lamp.

* Installation in tabletop position:

- Assemble the screws of the hinge as in Fig. 1.
- Assemble the wing nuts as in Fig. 1.
- Connect the plug into the mains.
- Press the switch to start the lamp.

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
Be sure that the voltage supply is the same as the one indicated on the characteristic plate of the equipment.

Do not use the equipment if it is NOT earthed.

If you have to change the plug, bear in mind the following:

Blue cable:	Neutral.
Brown cable:	Phase.
Yellow/green cable:	Earth.

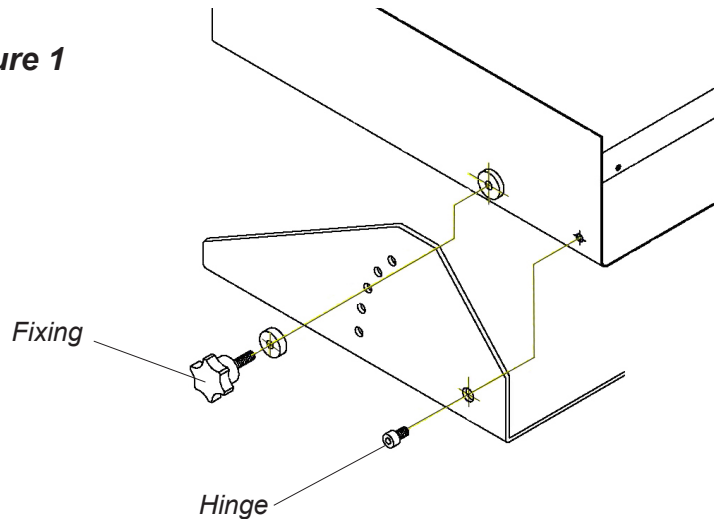Figure 1

GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the equipment by unathorized personnel by J.P. SELECTA, s.a.u. cancels the guarantee automatically.

MAINTENANCE

Before removing the cover, disconnect the equipment from the mains.

The manipulation of the internal electronic circuits of the equipment by unauthorized personnel can cause irreparable damage. Take it to one of the J.P. SELECTA's authorized technical services.

CLEANING:

For the cleaning of the different parts use the following products:

- Cleaning of stainless steel: Alcohol
- Cleaning of plastic: Alcohol with cotton duster.

SPARE PARTS

To guarantee the safety of the equipment, the spare parts must be purchased from J.P.SELECTA, s.a.u.

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