PLACA CIRCULAR DOBLE DUPLAC
DOUBLE CIRCULAR HOTPLATE DUPLAC

4000634
4000634

PLACAS CALEFACTORAS
HOTPLATES

INFORMACIÓN GENERAL

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the apparatus with liquids which can give off vapours capable of making explosive mixtures.*

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CONTENTS

GENERAL INFORMATION
CONTENTS
PACKING LISTS
EQUIPMENT DESCRIPTION
TECHNICAL FEATURES
INSTALLATION
OPERATION
MAINTENANCE
SPARE PARTS
GUARANTEE
ELECTRICAL DIAGRAM
"EC" CONFORMITY DECLARATION

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The standard equipment consist of the following components:

Pieza	Piece	Código / Code	Cant. / Qty.
Placa circular doble	Double circular hotplate	4000634	1
Manual de instrucciones	Instruction manuals	80111	1

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Double circular hotplate are designed to work with a big range of temperatures up to 400°C.

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The temperature regulation is done by independent energy regulator in each plate.

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Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

Código / Code	Nº de plazas / Nº of places	Ø Placa cm / Ø Plate cm	Medidas exteriores cm / Overall dimensions cm			Consumo por placa W / Consumption each plate W	Máxima temperatura °C / Maximum temperature °C	Peso Kg. / Weight Kg
			Alto / Height	Ancho / Length	Fondo / Width			
4000634	2	19	15	47	24	850	400	8.5

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Place the apparatus on a flat, horizontal, level surface, trying to leave a space of about 10 cm at the back and on each side.

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For a longer life of the apparatus provided with hotplate, it is recommended that its surface be covered as much as possible by the beaker to be heated. This avoid overheating and deformation of the plate by absorbing more the heat.

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Be sure that the voltage supply is the same as the one indicated on the characteristic plate of the apparatus.

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Do not use the apparatus if it is NOT earthed.

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If you have to change the plug bear in mind the following:

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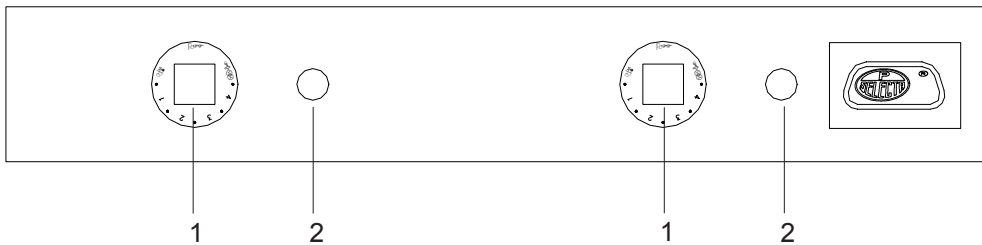
1. Move the control knob of temperature (1) from the "OFF" position (disconnected), to the "FULL" position (maximum power).
2. It will be noted that according to the control position, the ON-OFF cycle switch of the indicator lamp for the working of the heater (2) will be longer or shorter, which will permit adjusting to the desired working temperature.
3. To stop, move the control button to "OFF" position.

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WARNING!!! Caution with the hotplate surface, serious burns can be produced, even after switching off the apparatus.

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MAINTENANCE

Before removing the cover disconnect the apparatus from the mains.

The manipulation of the internal electronic circuits of the apparatus by unauthorized personnel can cause irreparable damage. Take it to one of the J.P.SELECTA authorized technical services.

CLEANING:

For the cleaning of the different parts use the following products:

Cleaning of stainless steel: Alcohol

Cleaning of plastic: Alcohol with cotton duster.

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GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use of causes beyond the control of J.P. SELECTA,S.A..

Any manipulation of the apparatus by unauthorized personnel by J.P. SELECTA,S.A. cancels the guarantee automatically.

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