ARMARIOS CONSERVADORES

UPRIGHT FREEZERS

TEMPLOW-S

2101278


INDICE

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INDEX

1 Safety	13
1.1 SAFETY ICONS	13
1.2 Risks that the operator can be subjected to	13
1.3 Qualified users	13
2 General information	14
3 Technical specifications.	14
4 Contents list	15
4.1 Standard contents	15
5 Introduction	15
6 Equipment description	15
6.1 Equipment	15
6.2 Features	15
6.3 Panel de mandos	16
7 Installation	16
8 Operation	17
8.1 Switching On the equipment.....	17
8.2 Brightness of the temperature display	17
8.3 Lock against accidental handling.....	17
8.4 Door alarm	18
8.5 Temperature alarm	18
8.6 Setting the temperature	18
8.7 Superfrost	18
8.8 Defrost	19
9 Maintenance	19
9.1 Cleaning	19
9.2 Changing over the door hinges	19
10 Guarantee	20

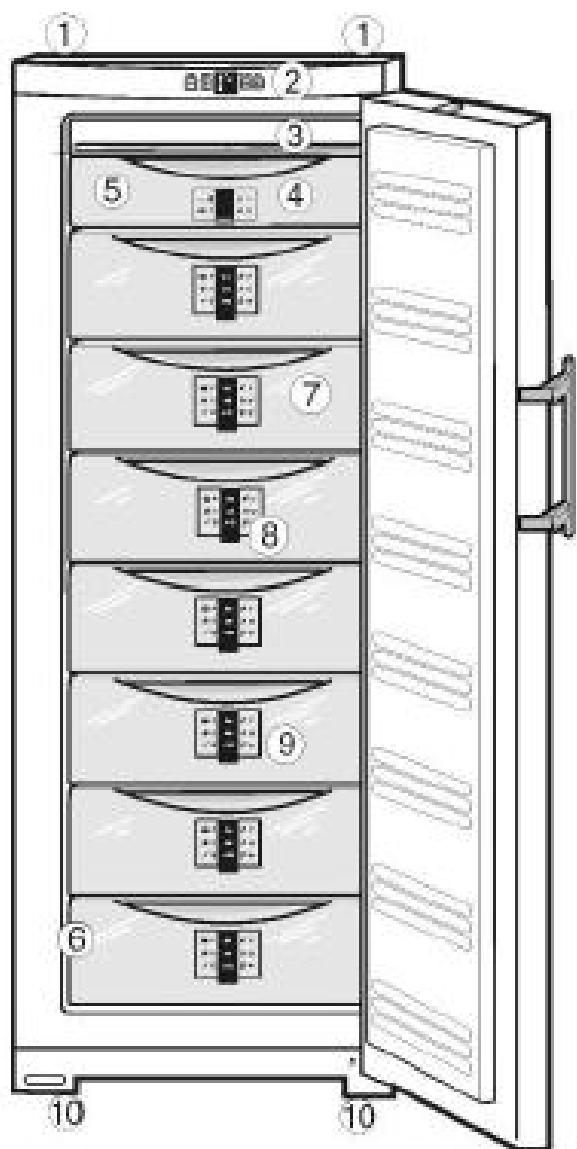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- 1) Handles for transport at the back.
- 2) Operating and control components.
- 3) NoFrost Device
- 4) Compartment
- 5) Freezing accumulator.
- 6) Nameplate.
- 7) Drawers.
- 8) Information system.
- 9) Vario Space
- 10) Adjusting feet, transport handles at the front, wheels at the rear.

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1 Safety

The equipment comes complete with safety features.

This manual indicates areas of possible risk.

1.1 SAFETY ICONS

Identifies risky situations and safety measures that should be adhered to.

The icons make reference within the paragraph, marked as a grey line.

Risk danger

Risk danger

Beware, follow instructions as indicated.

Electrical risk

Risk of electrical shock if accessing zones shown with this icon.

Beware, follow instructions as indicated.

Important information

- Important information in obtaining the best results or an optimum performance of the equipment.
- Important information to extend the life of the equipment and to maintain its optimum performance.

1.2 Risks that the operator can be subjected to

- Electrical risk.

1.3 Qualified users

This equipment should only be used by personnel that are qualified users.

This equipment should only be used by personnel that haS understood this manual or has previous instruction about the use of steam sterilizers.

2 General information

- 2.1 Handle the equipment with care. Unpack and check that all items coincide with the delivery note. If you see any discrepancy or damage of any kind, notify your distributor as soon as possible.
- 2.2 Read this manual before using the equipment.
- 2.3 These instructions are an important part of the equipment and should be made available to all users.
- 2.4 If in any doubt or clarity on the use of this instrument, contact your local distributor or the service department of J.P. Selecta, s.a.u.
- 2.5 **ATTENTION, AN EQUIPMENT NOT CLEAN AND DISINFECTED SHALL NOT BE REPAIRED.**
- 2.6 Any unauthorised modification, lack of maintenance or deliberate damage to the machine, transgresses the directive 89/655/CEE where the manufacturer will not be responsible for damages that could occur due to an undue care and attention.
- 2.7 Do not use the equipment for fluids that can produce vapours that can form inflammable or explosive mixtures.
- 2.8 Do not allow naked flames or ignition sources to enter the appliance.
- 2.9 Do not use electrical appliances inside (e.g. steam cleaners)
- 2.10 Avoid prolonged skin contact with cold surfaces or chilled/frozen food. This could cause pain, numbness and frostbite. Use protective gloves.

3 Technical specifications.

Voltage supply 115-230V 50/60Hz according to the characteristics plate indications.

Code:	2101278
Temperature range:	-15 to -32°C
Capacity:	257 litres
Dimensions:	
Height	184 cm
Width	60 cm
Depth	63 cm
Inner dimensions:	
Height	158 cm
Width	46 cm
Depth	43 cm
Power:	110 W
Weight:	74 kg

4 Contents list

4.1 Standard contents

The standard equipment consists of the following items:

- Templo-S 21012778
- Adjustement spanner
- Instruction manual 80115

5 Introduction

The upright freezer Templo-S is useful to conservation at low temperature. It has an external enclosure epoxy coated, an inner cabinet made from ABS chemical products-proof, changeable over door hinges, electronic digital temperature regulator, NTC probe, acoustic and luminous alarm of temperature, acoustic alarm disable key and superfrozen key.

6 Equipment description

6.1 Equipment


The upright freezer Templo-S is equipped with the following features:

- Digital temperature regulator
- NCT temperature probe
- 8 ABS drawers
- Acoustic and luminous alarm of temperature
- Acoustic alarm disable key
- Superfrozen key

6.2 Features

- Temperature regulation from -15°C to -32°C
- Changeable over door hinges
- Super frozen available
- Manual defrost with evaporating system

6.3 Panel de mandos

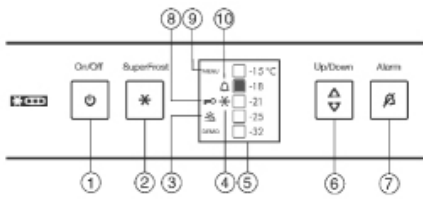1. ON / OFF key
2. SuperFrost key
3. Brightness Symbol
4. SuperFrost Symbol
5. Temperature indicator
6. Up / Down adjustment key
7. Alarm key
8. Block symbol for accidental handling
9. Menu symbol
10. Alarm symbol

7 Installation

7.1 Placement

Place the equipment near of a socket on a levelled floor (see «Power supply connection section) keeping in mind the following:

- When placing the equipment make sure that the pipes are not damaged.
- Once in position, use the provided spanner to adjust the feet so that the equipment is leveled.
- Do not place it in direct sunlight or near to a heating source.
- Do not place the equipment in wet places or where may get water splashes.
- The room where is installed the equipment has to be at a distance of, at least, 7 squared metres because, in case of refrigerant gas leakages, would not exist risk of flammable mixtures production.
- Before starting up, remove all the transport blocking.

DO NOT INSTALL THIS EQUIPMENT IN EXPLOSIVE ATMOSPHERES OR IN PLACES WHERE THERE ARE SOLVENTS OR OTHER PRODUCTS WHICH MAY PRODUCE EXPLOSIVE OR FLAMMABLE GAS.

7.2 Power Supply Connection

Make sure that the voltage given by the socket is the indicated on the feature label attached on the rear part.

The plug socket must be protected by a 10A to 16A fuse or circuit breaker.

Do not use the equipment unless it is correctly earthed.


If you replace the plug, keep in mind the following:

Blue wire:	Neutral
Brown wire:	Phase
Yellow/green wire:	Ground

8 Operation

8.1 Switching On the Equipment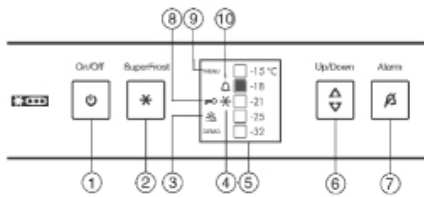

- Press the On/Off button(1).
- The appliance is switched on. The temperature display and Alarm symbol (10) flash until the temperature is cold enough.
- If "DEMO" is displayed, demo mode is activated.

8.2 Brightness of the Temperature Display

You can adjust the brightness of the temperature display to the light conditions of the room in which the appliance is installed.

8.2.1 Adjusting the brightness

The backlight can either be switched off or set to one of 5 levels. The backlight is switched off ex works.

- To activate the setting mode: Press the SuperFrost button (2) for about 5 seconds.
- The menu symbol (9) shines and the child lock symbol (8) flashes.
- Press the setting button (6) to select the brightness function.
- The child lock symbol (8) goes out and the brightness symbol (3) flashes.
- To confirm: Briefly press the SuperFrost button (2).
- The brightness symbol (3) shines.
- Using the setting button (6), select either "off" or the brightness wanted. The more display fields of the temperature display that light up, the brighter it is. If no field is lit up, it means "off".
- To confirm: Press the SuperFrost button (2).
- The brightness symbol (3) flashes.
- The brightness is adjusted to the new value.
- To deactivate the setting mode: Press the On/Off button (1) or wait for 5 minutes.
- The brightness symbol (3) and the menu symbol (9) go out.
- The temperature is indicated again in the temperature display.

8.3 Lock against Accidental Handling


The accidental handling lock function enables you to make sure that the equipment is not inadvertently switched off by an accidental handling.

8.3.1 To activate accidental handling lock

- To activate the setting mode: Press the SuperFrost button (2) for about 5 seconds.
- The menu symbol (9) shines and the accidental handling lock symbol (8) flashes.
- Briefly press the SuperFrost button (2) to select the accidental handling lock function.
- The accidental handling lock symbol (8) shines. The -15°C and -21°C LEDs shine in the temperature display.
- Briefly press the SuperFrost button (2) to activate the accidental handling lock.
- The accidental handling lock symbol (8) flashes. The -15°C and -21°C LEDs go out.
- To deactivate the setting mode: Press the On/Off button (1) or wait for 5 minutes.
- The menu symbol (9) goes out and the temperature is displayed again in the temperature display. The accidental handling lock symbol (8) shines.

8.3.2 To deactivate accidental handling lock

- To activate the setting mode: Press the SuperFrost button (2) for about 5 seconds.
- The menu symbol (9) shines and the accidental handling lock symbol (8) flashes.
- Briefly press the SuperFrost button (2) to select the accidental handling lock function.
- The accidental handling lock symbol (8) shines. The -18°C LED shines in the temperature display.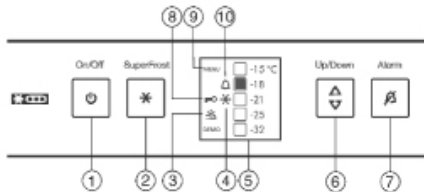

- Briefly press the SuperFrost button (2) to deactivate the accidental handling lock.
- The accidental handling lock symbol (8) flashes.
- To deactivate the setting mode: Press the On/Off button (1) or wait for 5 minutes.
- The menu symbol (9) goes out and the temperature is displayed again in the temperature display. The accidental handling lock symbol (8) no longer shines.

8.4 Door Alarm

If the door is open for longer than 60 seconds, the audible warning will sound. The audible alarm is automatically silenced when the door is closed.

8.4.1 Muting the door alarm

The audible alarm can be muted when the door is open. The sound switch-off function is active as long as the door is left open.

- Press alarm button (7).

8.5 Temperature Alarm

The audible alarm sounds if the freezer temperature is not cold enough. The temperature display and the alarm symbol (10) flash at the same time. The cause of the temperature being too high may be:

- warm fresh product was placed inside
- too much warm ambient air flowed in when rearranging and removing product
- power failure for some time
- the appliance is faulty

The audible alarm is automatically silenced, the alarm symbol (10) goes out and the temperature display stops flashing when the temperature is sufficiently cold again.

If the alarm status persists: (see Malfunction).

Note:

products may be spoilt if the temperature is not cold enough.

- Check the quality of the product.

8.5.1 Muting the temperature alarm

The audible alarm can be muted. When the temperature is sufficiently cold again, the alarm function is active again.

- Press alarm button (7).
- The audible alarm is silenced.

8.6 Setting the Temperature

Temperature setting to be recommended: -18°C.

The temperature can be changed continuously. Once the -32°C setting is reached, it starts again with -15 °C.

- To access temperature adjustment: press setting button (6) once.
- The LED of the current temperature flashes in the temperature display.
- Press the setting button (6) the number of times needed until the required temperature shines in the LED display.


Note:

- Long pressing of the setting button sets a slightly colder value within a small temperature range (e.g.: between -15°C and -18°C). The LED of the next lower temperature range then shines in the temperature display.

8.7 Superfrost

This function serves to freeze the products more quickly. The maximum freezing capacity of the Templo-S is 20kg of fresh products at room temperature for every 24 hours.

- Briefly press the SuperFrost button (2).
- The SuperFrost symbol (4) shines.
- The freezer temperature drops, the appliance operates with the maximum refrigerating capacity.

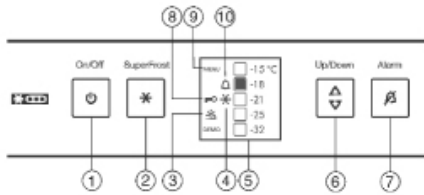In case of a small amount of product to be frozen:

- wait approx. 6 h.
- Place wrapped produce in the deep bottom drawers.

In case of the maximum amount of product to be frozen (see type plate):

- wait about 24 h.
- Remove the deep bottom drawers and place produce directly on the bottom shelves.

- SuperFrost is automatically deactivated after about 65 h.
- The SuperFrost button (4) extinguishes once freezing is completed.
- Place product in the drawers and push the latter back in again.
- The appliance continues to operate in the energy-saving, normal mode.

8.8 Defrost

The No-Frost system is responsible for the automatic defrosting of the appliance.

Moisture from the freezing process adheres to the evaporator. It will be defrosted and periodically evaporated automatically.

Note:

- Do not defrost manually.

9 Maintenance

9.1 Cleaning

- BEFORE START TO CLEAN UNPLUG THE EQUIPMENT FROM THE SOCKET
- Clean the cabinet with warm water mixed with a small quantity of the soap.
- DO NOT USE CLEANING STEAM DEVICES. IT IS A DANGER FOR THE PEOPLE AND FOR THE EQUIPMENT.

- Try the water used do not reach the air inled grid or the electric elements.
- Dry the equipment with a soft dry cotton cloth.
- Clean once a year the duster accumulated on the compresor and the rear part grid. This acumulation increases the energy consumption.
- During the cleanig consumption, pay attention on avoiding that pipes or other parts may get broken.
- After finishig the cleaning, connect the freezer again, activate the function superfrost and put inside the low freezed products as soon as the temperature reaches the set temperature again.

If the appliance is to be left out of operation any length time, empty it, disconnect from the mains and clean it. If necessary, left the door opened in order to avoid unpleasant smells.

9.2 Changing over the Door Hinges


To change the opening direction of the door, do as follows: disconnect the equipment from the mains and ensure that the following tools are to hand:

- Torx 25
- Torx 15
- Open-ended spanner size 6
- Screwdriver
- Open-ended spanner included in supply package

ATENCIÓN

Risk of injury if the door tips!

Take good hold of the door and set down the door carefully.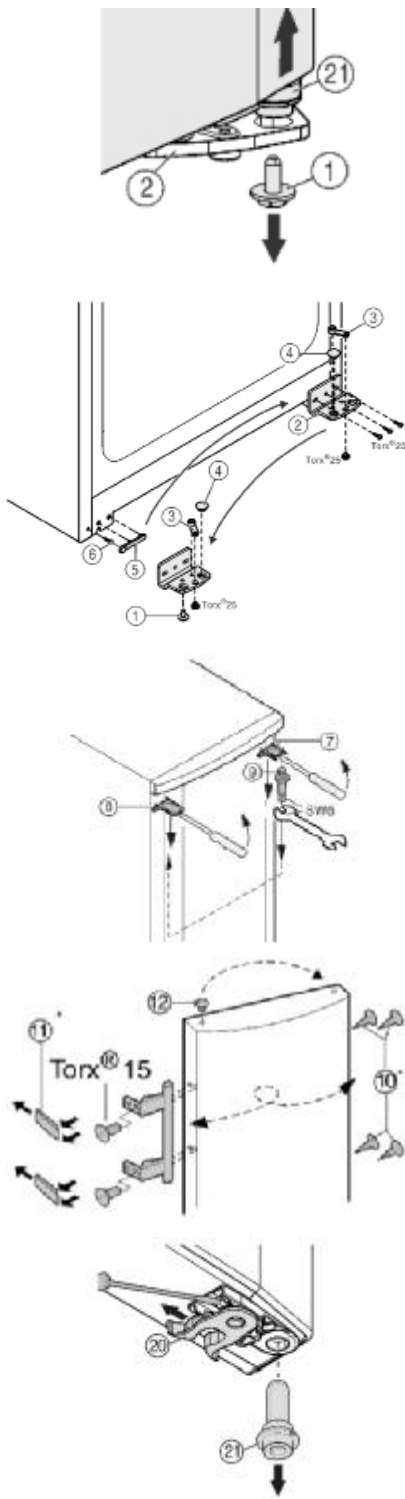

- Unscrew the securing screw on the bottom right turn hinge.
- Open the door.
- Hold the door on the handle side and at the bottom and lift it.
- The bearing pin (21) should come out of the turn hinge (2).
- If the bearing pin (21) does not come out, press the pin out from underneath.
- Swing door out at the bottom and remove it.
- Unscrew the turn hinge (2).
- Unscrew the bearing element, (3) transfer it to the opposite location hole of the turn hinge and screw it firmly into place.
- Take the stopper (4) out of the turn hinge and transfer it to the receiving hole on the opposite side of the turn hinge.
- Carefully lift off the cover (5) on the handle side.
- Undo the screw (6) and transfer it to the opposite side.
- Put the cover (5) back on.
- Screw the turn hinge (2) firmly into place on the new hinge side (with 4 Nm), possibly using a power screwdriver.
- At the top, use a screwdriver to unclip cover (7) and cover (8) at the front, and remove by pulling out and down.
- Unscrew the bearing pin (9) and screw it firmly into place on the opposite side (with 4 Nm).
- Refit the cover (7) on the side of the bearing pin: insert at the back, snap in at the front.
- Fit the cover (8) on the opposite side: insert at the back, snap in at the front.
- Lift the stopper (12) out of the door bearing bush and transfer it.
- Detach door handle, stoppers (10) and pressure plates (11) and transfer to the opposite side.
- When mounting the pressure plates, make sure they snap into place properly.
- Transfer the spring clamp (20): Depress the latch nose and pull the spring clamp off over it.
- Slide the spring clamp into place on the new hinge side until it clicks into place.
- Pull the bearing pin (21) out of the door bush and transfer it to the opposite side together with the washer. The latch nose must point to the inner side of the door, the notch to the outer side.
- Hang the door on the bearing pin (9) at the top.
- Swing the door in at the bottom and insert the bearing pin (21) into the turn hinge. If necessary, turn the bearing pin until it snaps into place.
- Insert the securing screw (1) at the bottom into the bearing pin and tighten it (with 4 Nm).
- If necessary, align the door flush with the appliance housing using the two slots in the turn hinge. To do so, unscrew the middle screw.

WARNING

Risk of injury due to the door dropping out!

If the bearing parts are not screwed into place firmly enough, the door may drop out. This may lead to severe injuries. What is more, the door may not close and therefore the appliance may fail to cool properly.

- Screw on the bearing brackets/bearing pins tightly (4 Nm).
- Check all of the screws and retighten if necessary.

10 Guarantee

This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the equipment by personnel not authorized by J.P. SELECTA, s.a.u. automatically cancels the guarantee.