EF-1425

4001425

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0. GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact your supplier or J.P. Selecta's technical service.*
- 5) ***IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.***
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*

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2. PACKING LIST

The standard equipment consists of the following components:

Componentes	Components	Código/Code	Cant. / Qty.
Mueble soporte unidad	Support furniture unit	9457	1
Rampa	Ramp glass	47030	2
Colector	Collector glass	47031	1
Frasco PYREX 1L SVL 42	Flask PYREX 1L SVL 42	47033	1
Tapones completos SVL 42	Stoppers SVL 42	28323	2
Capuchón unión SVL 42	Union SVL 42	28310	7
Capuchón unión SVL 22	Union SVL 22	28263	2
Anillo estanqueidad 30Ø int	Sealing ring 30Ø int	21168	1
Anillo estanqueidad 14Ø int	Sealing ring 14Ø int	30149	2
Juntas tóricas 34x4 viton	Gasket 34x4	21190	6
Manguera vitón 12 Ø(negra)	«Viton» tube 12 Ø(black)	46085	1 m
Manual de instrucciones	Instructions manual	80145	1

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6. INSTALACIÓN

4. FEATURES

Stainless steel support AISI 304.
Ramp and carries crucibles in discussed glass.
Modular construction that facilitates its cleanliness.

5. DESCRIPTION

Specially designed unit to accomplish the cold extraction with acetone or other solvents, of 6 samples simultaneous.

It is formed by a container of 1litro (in the one which be recovered the solvent) and 6 adapting to fix the filtering crucibles .

Need a water jet pump connected to the suction pipe to accomplish the extraction.

To work with a number samples minor is necessary put complete stoppers (supplied with the unit) in the empty localities.

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Put aprox. 60mm of VITON tube (black) in the ramps ends.

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Put the union SVL22 and the sealing ring 14 Ø int.

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Put the ramps on the unit support furniture.

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Put the union SVL42 and the sealing ring 30 Ø in the neck of the collector.

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Put the collector on the flask PYREX 1L SVL42.

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Locate the stainless steel support in a stable and horizontal flat surface.

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Put the free ends of the VITON tube (black) in the upper orifices of the collector.

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Connect a "viton" tube (black) end to the water jet pump and install it.

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Connect the other end of the tube of the water jet pump to the collector.

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Put the gasket 34x4 and the union SVL42 on the crucibles that we want to use.

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Put the crucibles on the ramps and the stoppers SVL42 on the orifices without sample. Now the unit is ready to be used.

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7. INITIAL CHECKING

Add 25 ml of water in each filtering crucibles and open the water current through the water jet pump. Check that the filtration is produced quickly. Otherwise, check that all the stoppers and unions are well tightened.

To obtain a most quick filtering, increase the water circulation through the water jet pump.

8. MAINTENANCE

It is advisable to accomplish an exhaustive cleanliness after each use.

Dismantle the components of the unit and insert them in a ultrasonic bath with distilled water and an appropriate detergent (BIO-SEL) for around 10 minutes.

Rinse with abundant distilled water and let it dry.

Before each use, make an inspection of the tubes and replace them if they are damaged.