DESTILADOR DE AGUA PARA AUTOCLAVES
WATER DISTILLER FOR AUTOCLAVES

4001729

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1 Packing list

The distiller is supplied with the following items:

- | | |
|-------------------------|-------------------|
| 1. Distilled water tank | 2. Water guiding. |
| 3. Filter (2 units) | 4. Power cord. |

2 Installation

Description:

Benchtop water distiller for special usage in autoclaves.

Suitable location:

- Place on a flat level and stable surface, suitable for the unit weight.
- Do not cover the ventilation holes.
- Place it near to power socket.

Connection to power supply:

- Check that voltage on socked match the one indicated on the plate.
- Do not manipulate the power cords with wet hands.
- Power socked must be provided with protective ground terminal.

3 Technical specification

Power consumption:	750W
Distillation rate:	2 Litres/h
Vessel volume:	4 Litres
Power supply Voltage:	230V AC \pm 10% (50/60Hz) See plate
Weight:	3.6Kg
Dimensions:	245 x 285 x 512 mm

4.1

4.2

4.3

4 Operation

- Unplug the header. Fig 4.1 and remove it. Fig 4.2
- Fill the internal vessel, with water up to «FULL». Fig 4.3
- Check that filter is properly fixed.
- Fix the header. Fig 4.2 and plug the wire. Fig 4.1
- The fan starts to run to cold the condenser.

Note: If the fan does NOT start, press «SPSP».

- When reaching the boiling temperature, the distilled water starts to flow through the faucet.
- When the water is finished, the distiller stops automatically.
(It is suggested to unplug from the mains when the unit is not used)
- Wait near 30 minutes to let the unit become cold, before removing the header and clean the internal vessel and header bottom.

Tip: Distillation time can be shorten if hot water is used.

5 Cleaning

The water tap distillation left calcium residues on the vessel walls. This residues affect the distillation rate and could let the unit out of order if they are not regularly cleaned.

- Use an anti- calcium, applied with a non-abrasive cloth to clean the inner parts.
- Clean each 2 - 3 distillations. Do not let accumulate the calcium residues on the inner surfaces.
- Deionized water can be used to avoid calcium residues.
- Change the output filter every 3 moths. (Assuming a daily usage)

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Electrical equipment for measurement, control and laboratory use. EMC Requirements.

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Safety requirements for electrical equipment for measurement, control and laboratory use. Part 1: General requirements.

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7 Warranty

The warranty covers the defective components or faulty assembling of the unit for 1 year from the purchasing date. Keep purchasing bill as the only valid document for the purchasing date.

The warranty does not covers the damages caused for a unit malfunction.

Return goods are not accepted without previous J.P. SELECTA,s.a. approval.