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1 Introduction

Ultrasonic cleaning baths are usually used in:

General laboratory for:

- chemical reactions acceleration
- liquids degassing
- cell disruption
- sieves cleaning, pipettes, micro-pipettes, cuvettes, viscosimeters
- decomposition of radioactive substances.

Odontology for:

- prosthesis and instruments cleaning.

Optics: for frames, lenses and contact lenses cleaning.

Medicine: for instruments in general cleaning, forceps, probes, scalpels.

Jewellery and watch making: for any kind of watches and jewels cleaning.

Industry: for all electronic components cleaning, printed circuits board, semiconductors, petrol injectors, sieves and filters.

2 Working Principle

Ultrasonic cleaning consists of the use of high frequency sound waves (40 KHz) which produces millions of low pressure microscopical bubbles with high penetration, eliminating impurities, polluting agents and dirt from the parts or material which must be cleaned.

3 Technical Specifications

Model	Inner dimensions (HxWxD) mm	Capacity	Generator power	Heating power	Temperature range	Adjustable time	Drainage
		L	W	W	°C	min	
3000865	150 x 240 x 140	5	60	200	0-80	1-99	Yes
3000866	150 x 300 x 155	6	70	400	0-80	1-99	Yes
3000867	150 x 300 x 240	10	150	400	0-80	1-99	Yes
3000868	150 x 500 x 300	22	450	900	0-80	1-99	Yes
3000864	200 x 500 x 300	30	800	900	0-80	1-99	Yes

4 Packing List and Accessories

Packing List:

Ultrasonic bath
Wire basket
Lid with handle
Drain plug
Power cord
Instructions manual

Accessories (not included on the standard package):

Bio-Sel cleaning solution: Code 1000025

5 Warnings

1. Read carefully this manual before using the machine.
2. The grounding device must be connected before using the equipment.
3. Do not use the bath with flammable liquids or those that could give off vapours capable of making explosive mixtures.
4. Take care of using a cover when using volatile solutions.
5. Don't let the temperature of cleaning solution reach 80°C.
6. Never start the machine when there's neither water nor cleaning solution.
7. This appliance is intended for professional use in the laboratory.

6 Operation

1. Power socked must be provided with protective ground terminal.
2. Fill bath with water, at least 3/4 full.
Use demineralized water.
Use a suitable detergent according to the material to be cleaned.
3. Switch the equipment on.
4. Set the temperature, according to the cleaning requirements.
5. Place articles to be washed into the basket and set the cleaning time. Normally, time can be set between 3 and 30 minutes, and can be prolonged for tools which are especially difficult to clean.
6. For heavy articles, suspend them through a hanger in the cleaning solution to increase the cleaning effect and the cleaning quality.
7. If you don't use the equipment for a long period of time, do not leave the tank full of liquid.
8. Empty the cleaning solution and clean the bath tank.

7 Temperature and Time Selection

1. Set the ultrasonic cleaning time to max. 99 minutes. Increase or reduce the time by means of the keys till you set the needed cleaning time.
2. Adjust the temperature (0-80°C). Increase or reduce the temperature by means of the keys until you set the needed temperature for the cleaning solution.
(Note: the bath only heats, it is not able to cool)

8 Service and Maintenance

This equipment should only be manipulated by authorized maintenance personnel.

A periodical checking should be done to the equipment.

1. Visually check if there is water leakage or dampness.
2. Visually check if the power connection cable is in good condition.
3. Visually check if the cooling rear fan accumulates dust on it. Clean with a brush or with a vacuum cleaner.

9 Warranty

This product is guaranteed for one year since the purchasing date. The guarantee does not cover damage caused by incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the equipment by unauthorized personnel by J.P. SELECTA, s.a.u. automatically cancels the guarantee.

Notice to customers:

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.

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**ATTENTION:****DO NOT START UP THE BATH WITHOUT WATER ON THE TANK.**

When switching on the main switch («ON/OFF»), the heater will be activated.

The tank heater is independent of the «START/STOP» key and it can be switched on directly by the main switch.

If heater is activated without water on the bath tank, the equipment could be damaged.