ELEVADOR PARA AUTOCLAVE VERTICAL

CRANE FOR VERTICAL AUTOCLAVE

1002426

1. General Information

Handle the parcel with care and check that contents coincide with the packing-list. If any part is damaged or missing, please advise the distributors immediately.

Do not install or use the equipment without reading this handbook first. This instructions manual must be attached to the equipment and must be available for all users.

If you have any doubts or enquiries, contact with J.P. SELECTA's technical service.

Do not use the equipment in explosive, flammable or corrosive atmospheres.

2. Packing list & accessories

The crane is supplied together with 6 screws for its attachment to the autoclave.

3. Technical specification

Reference code:	1002426
Dimensions: maximum high in up position	2400 mm
Volage supply:	230 VAC 50Hz
Power consumption:	70 W (0.5A)
Weight:	23 Kg
Maximum load:	50 Kg

4. Information about safety in the use of the equipment

4.1 Staff qualification

This equipment may only be used by properly trained personnel who have read and understand these instructions or have been properly trained in this equipment operation.

4.2 Use with hazardous substances

Do not treat flammable or explosive materials or those which chemically react with violence.

4.3. Maintenance

Unplug the mains before removing covers to handle inside.

The manipulation of the equipment internal electronic circuits by unauthorized personnel can cause irreparable damage. Contact with J.P.SELECTA's authorized service.

Use glass cleaner or similar to clean the accessible parts. Do not use aggressive solvents.

4.4. Warranty

This product is guaranteed for one year. It does not cover damages caused by an incorrect use or causes beyond the control of J. P. SELECTA, s.a.u.

Any manipulation of the equipment by personnel not authorized by J.P. SELECTA, s.a.u. automatically cancels the guarantee.

5. Equipment description

Crane 1002426 has been designed for loading and unloading J.P. SELECTA's autoclaves baskets.

It is suggested to use a transport trolley between 70 and 80 cm high to transport the baskets up to the crane reach.

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6. Intallation

The crane 1002426 only could be assembled on the autoclaves provided with the fixing holes at the left hand side.

A power plug will be needed near the crane.

Fix it with 6 screws and nuts according the following drawing.

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This part could be reversed to operate with low profile baskets.

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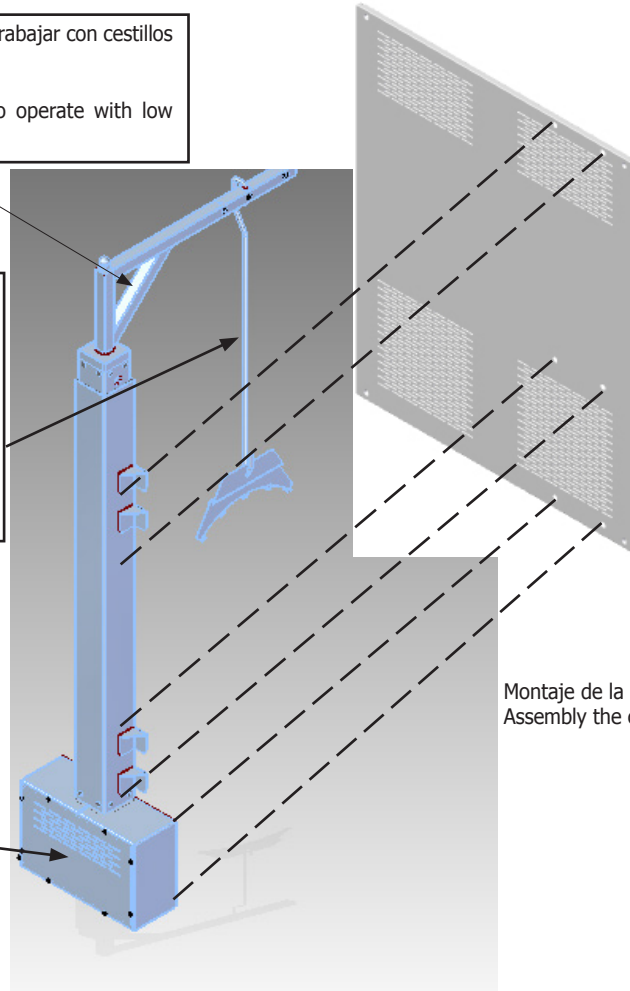
Guide with hand to avoid hits on the basket.

Al quitar la tapa.
If cover is removed.

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For crane operation, we suggest to guide the basket hang with one hand and to operate the up/down switch with the other.

Use a transporter from 70 to 80cm high to load /unload the baskets.

If baskets have a low height, the top hang support could be assembled at reverse position. (See figure 1)

J.P. SELECTA, s.a.u. 2019