2000963	36L
2000964	80L
2000965	150L

Prebatem-TFT

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1. General Information

- Handle the package with caution. Unpack and check that the contents match what is indicated in the "Packing List" section. If a damaged component or the absence of any is observed, please advise the distributor immediately.
- Do not install or use the equipment without reading this instruction manual before. These instructions are an inseparable part of the equipment and should be available to all users.
- Check any doubt with J.P. SELECTA's technical service.
- Do not use equipments in explosive, flammable or corrosive atmospheres.

2. Packing List

Model:	2000963	2000964	2000965
Inner volume:	36L	80L	150L
Shelves:	2000022 x2	2000023 x2	2000025 x2
Guides:	2000012 x4	2000013 x4	2000015 x4

3. Technical Specifications

Model:	2000963	2000964	2000965
Inner volume:	36L	80L	150L
Inner dimensions: height/width/depth (mm)	400/300/300	500/400/400	500/600/500
External dimensions: height/width/depth (mm)	600/650/490	700/750/590	700/950/680
Power voltage supply:	230 VAC 50Hz		
Electrical power (W):	310		
Weight: (Kg)	54	73	94
External fuse:	Replaceable (Type: 10x5mm)		
Temperature range:	5.0°C ... 60.0°C		
Display resolution:	0.1°C		
Temporized operation:	Weekly, hour, minute		
Temperature profiles:	Up to 10 profiles of 6 segments. Segment from 1m to 99 hours		

3.1 Communications

Data logging:	USB pen drive
ETHERNET:	Control panel remote access
Serial:	RS-232
Wi-Fi:	Optional
Bluetooth:	Optional

CAPABILITIES	SPECIFICATIONS	
	at 10°C	at 37°C
Stability	±0.1°C	±0.1°C
Homogeneity	±0.3°C	±0.2°C
Set error	±0.4°C	±0.2°C

4. Information about safety in the use of the equipment

Hazard risks

Electrical risks

Risk of high temperature contact

4.1 Risks the operator is subjected to

- Risk of burns
- Electrical risks

4.2 Staff qualification

This equipment may only be used by properly trained personnel who have read and understood these instructions or have been properly trained in this equipment operation.

4.3 Use with hazardous substances

Do not treat with flammable or explosive materials or with those which chemically react with violence.

4.4 Maintenance

Unplug the equipment from the mains before removing the oven's cover to handle inside.

The manipulation of the equipment internal electronic circuits by unauthorized personnel can cause irreparable damage. Take it to one of J.P. SELECTA's authorized technical services.

It is recommended to clean the accessible parts of the oven with a cloth and a cleaner, or a similar solution. Do not use aggressive solvents.

4.5 Warranty

This product is guaranteed for one year. It does not cover damages caused by an incorrect use or causes beyond the control of J. P. SELECTA, s.a.u.

Any manipulation of the equipment by personnel not authorized by J.P. SELECTA, s.a.u. automatically cancels the guarantee.

5. Equipment Description

PREBATEM ovens are used in sample thermal treatments, for its desiccation, cured, culture or conservation at a constant temperature. The working range includes temperature at 20°C which is very useful to temper and prepare measurement and calibration standards.

Heating and cooling system by semiconductors with electronic regulation and digital reading. The cooling system is affected by ambient temperature. The minimum temperature increases with the ambient temperature according to the following table:

Ambient temperature	Minimum temperature
25°C	5°C
26°C	6°C
27°C	7°C
28°C	8°C
29°C	9°C
30°C	10°C


The oven allows programming temperature profiles to optimize heat treatment.

The equipment supports several communication types for data logging and remote control to monitor the process.

1. Main TFT screen
2. Main switch
3. Safety thermostat
4. USB connection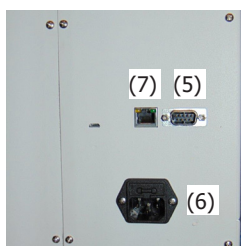

5. RS232 connection
6. Mains power connector + fuse
7. ETHERNET connection

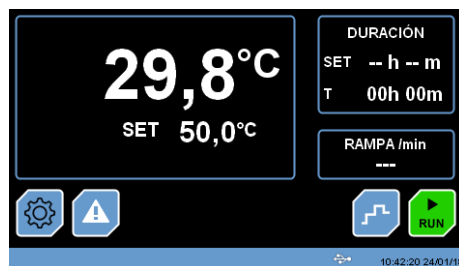
Fig. 6.3.1 MAIN SCREEN (START)

6. Installation

Once the equipment is unpacked, visually check that it is in good conditions. Keep the original packing for a few days.

6.1 Location

Place the machine on a flat level and stable surface, suitable for the weight of the equipment.

For comfort, leave a free space of 50cm around the equipment and a minimum of 20cm at the back.

6.2 Connection to the mains

Choose a close electrical outlet and appropriate for the equipment power. Use the supplied connection cable or a similar one.

For your safety, the socket must be grounded.

Before plugging, check that the voltage and the frequency of the electrical supply corresponds to the one indicated on the label of the equipment.

6.3 Starting

Choose an electrical outlet near the equipment and appropriate for the equipment power. Use the supplied connection cable or a similar one.

The oven starts with the O/I switch (2).

It starts with the **MAIN SCREEN**. (Fig. 6.3.1.)

(**RUN** & **STOP** buttons (icons) are activated according to the state they were when switching the oven off the last time with the O/I switch.)

From this screen, you can access the different operating options:

Settings menu. See **8 Settings**.

Simple work mode / programming. See **7 Operation**


Start / stop the oven.

Status bar. It indicates in which menu it is located, the date and time and if there is any type of communication connected (in the example, the USB connection icon appears).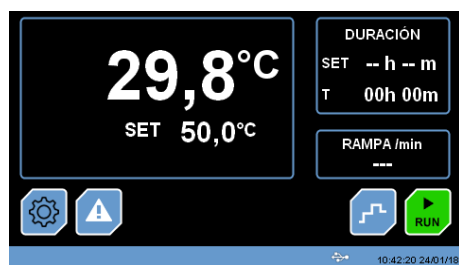

MAIN SCREEN (START)

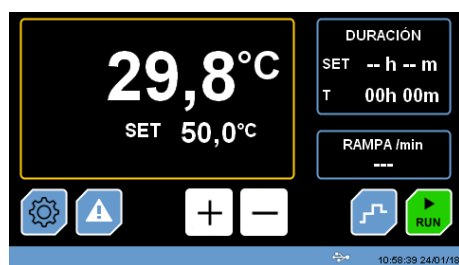
Fig 7.1 Screen to set temperature and time.

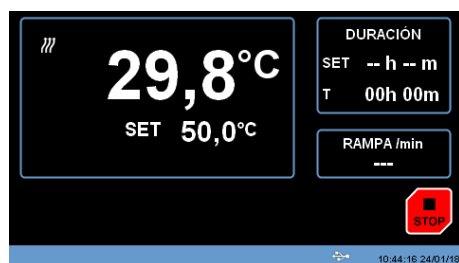
Fig 7.2 Screen to stop a cycle.

7. Operation

All the equipment functions are established through the touch screen. Press gently with your finger in the center of each icon.

This screen shows the status of the oven before disconnecting it by means of the main switch (0/I).

This allows the oven to resume its task in the event of a power cut.

7.1 Set temperature and working time

For most applications, this is the usual work method. If you need to create temperature profiles or program the start-up of the oven automatically, follow the steps in section 7.3.

To select the working temperature:

- From the **MAIN SCREEN**, click on the value to be modified. The frame of the modifiable parameter will be boxed in yellow. In the example, the temperature can be modified.

The **+** and **-** keys appear to increase / decrease the value of the adjustment.

- After a few seconds without pressing any key, the yellow frame disappears and the adjusted value is selected.

- The same operation can be performed to modify the working time or the temperature ramp.

A time of **--h --m** establishes a continuous operation.

7.2 Start / Stop the cycle


- From the **MAIN SCREEN** press:

RUN to start the cycle (the icon will change its appearance with the STOP message).

STOP to stop the cycle (the icon will change its appearance with the RUN message).

Note: When the duty cycle is in operation, a heat or cold symbol appears depending on the type of energy that the heating element is generating.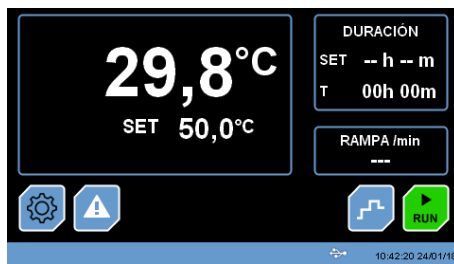

MAIN SCREEN (START)

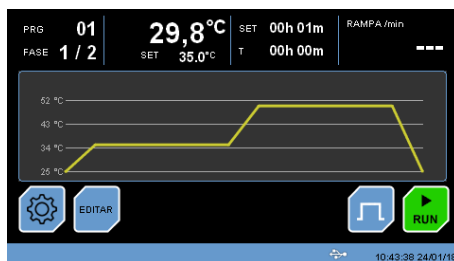
Fig. 7.3 Temperature profiles

Fig. 7.4 Editing temperature profiles

Fig. 7.4.1 Modify temperature, ramp and time parameters.

Fig. 7.4.2 Modify maximum and minimum temperature limits.

7.3 Programming temperature profiles

A temperature profile is a sequence of steps defined by a temperature and a time. It also allows you to define a schedule for the start / end of the process.

The equipment allows defining up to 10 profiles of 6 steps each.

From the main screen, press this button to work in "Programming" mode; the temperature profile will appear graphically as shown in Fig 7.3.

To start the program, simply press the **RUN** key (the button will change its appearance with the **STOP** message).

If during the program operation you wish to stop the process, press the **STOP** button.

From screen 7.3, click on this button to return to the main screen.

7.4 Editing temperature profiles

From screen 7.3, press the **EDIT** button to select and modify the program. The screen will show the appearance of Fig. 7.4.

Press the **EDIT** button to modify the temperature, time and ramp parameters, as shown in Fig 7.4.1.

Press the **ALARMS** button to define the ranges between which the system should work. If the reading of the temperature is above or below the defined values, a warning message will appear telling that the programmed limits have been exceeded. These alarms are only informative, they do not stop the operation of the equipment.

Press the **CYCLICAL** button to repeat indefinitely the selected program.

7.4.1 Set temperature, time and ramp parameters

The **+** and **-** keys allow you to add steps to the profile and the left / right arrows allow you to move between the different phases of the profile.

Press the **BACK** button to return to the main editing screen of profile 7.4.

7.4.2 Set maximum and minimum temperature limit values

If you want the oven to show an alarm message when the temperature is outside margins, set a value for over temperature and for low temperature.

Press on the value and use the **Up/Down** keys to adjust these limits.

Press the **BACK** key to return to the main editing screen of profile 7.4.

Fig 7.4.3 Selection date and time of program start


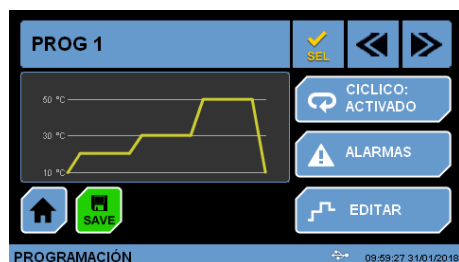
Fig. 7.4.4 Selection of working program

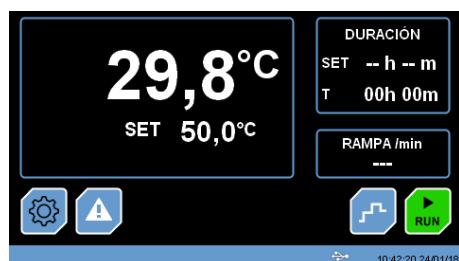MAIN SCREEN (START)

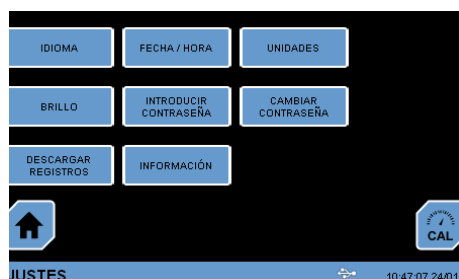
Fig. 8 SETTINGS SCREEN

7.4.3 Selection of date and time of program start

From screen 7.4.3 you can plan the week days and the start time of the cycle.

Press "Activate" to activate programming.

Press the **BACK** button to return to the main editing screen of profile 7.4.

NOTE: If a program is running, the cycle will not turn off. It has to be stopped.

7.4.4 Selection and saving working program

To select one of the 10 programs, press the **SEL** button from screen Fig.7.4 to select the working program of the oven.

Press the **SAVE** button to save all the parameter changes established in the program.

8 Settings

Press this button from the main screen to enter the settings screen as shown in Fig 8.

Modifying the settings parameters can affect the equipment operation. Read carefully the entire chapter so that the equipment works according to the process needs.

Most sensitive temperature control parameters are protected by a password. See the corresponding chapter.

Parameters settings are grouped according to their function as shown below:

LANGUAGE

Allows you to select the screen language.

DATE/TIME

Adjusts date and time of the oven. This adjustment is important if you work with programming.

UNITS

Sets parameters units.

BRIGHTNESS

Sets brightness of the screen backlight.

ENTER PASSWORD

Enters the password for accessing the calibration menu.

CHANGE PASSWORD

Changes initial factory password.

ALARMS

Sets margins out of which the oven gives an error.

LOGS DOWNLOAD

Press this button and the control will allow downloading of data stored in memory to a USB memory connected in the front part.

INFORMATION

Shows the firmware version and gives access to the menu for its update.

Fig. 8.1 Language selection

8.1 Language

When entering this setting, screen Fig 8.1 appears. Use the **Up/Down** arrows to select the language you want to appear in all menus.

Press the **BACK** key to return to the settings menu or the **HOME** key to return to the initial screen.

Fig. 8.2 Date / Time selection

8.2 Date/Time

Screen Fig 8.2 appears when you enter this setting. Click on the parameter to be modified and use the **Up/Down** arrows to select the corresponding value.

Press the **BACK** key to return to the settings menu or the **HOME** key to return to the initial screen.

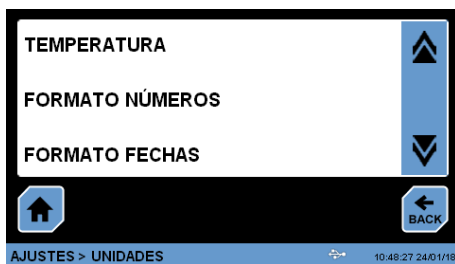
Fig. 8.3 Units selection

8.3 Units

When entering this setting, screen Fig 8.3 appears. Use the **Up/Down** arrows to select the parameter from which you want to change the unit.

Fig. 8.3.1 Temperature unit selection

Fig 8.3.1 shows the screen for selecting the temperature units. By using the **Up/Down** arrows you select between °C and °F. If this value is changed, all displayed values will be referenced to this unit.

Press the **BACK** key to return to the settings menu or the **HOME** key to return to the initial screen.

Fig. 8.4 Setting the screen brightness

8.4 Brightness

When entering this setting, screen Fig 8.4 appears. Use the **Up/Down** arrows to select the brightness percentage of the screen.

Press the **BACK** key to return to the settings menu or the **HOME** key to return to the initial screen.

Fig. 8.5 Setting alarm margins

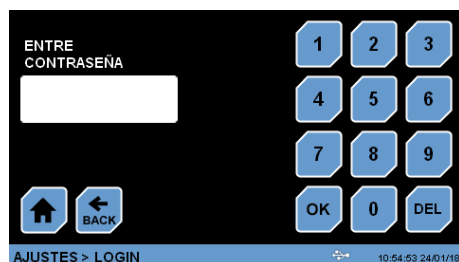
Fig. 8.6 Password entry

Fig. 8.6.1 Temperature Offset modification

Fig. 8.7 Password modification

Fig. 8.8 Logs download

8.5 Alarms

When entering this setting, screen Fig 8.5 appears. Use the **Up/Down** arrows to select the value of over temperature or low temperature you want for the temperature range.

This setting is simply a warning, it does not stop the operating cycle. This notice will be reflected in the memory log.

Press the **BACK** key to return to the settings menu or the **HOME** key to return to the initial screen.

8.6 Temperature Offset Adjustment

The Offset adjustment is protected by a password to prevent any user from modifying this parameter.

Press the **CAL** button to access this setting.

If the **CAL** button is pressed without entering the password, it will indicate that a valid password value is necessary to continue with the action. Therefore, it is necessary to select **ENTER PASSWORD** option before pressing the **CAL** button.

The default password value for the calibrator is "0000".

Press on the temperature Offset parameter and use the **Up/Down** keys to modify this value.

Press the **BACK** key to return to the settings menu or the **HOME** key to return to the initial screen.

The password value remains valid until the oven is switched off and on again. Perform this action to prevent another user from modifying the value.

8.7 Password Change

If you want to change the factory password, use this menu and follow the instructions that appear to change the password.

Press the **BACK** key to return to the settings menu or the **HOME** key to return to the initial screen.

8.8 Logs Download

To download the data stored in the control, select the option **LOGS DOWNLOAD**. Screen Fig. 8.7 will appear.

Connect a memory pen drive to the front connector (4) and follow the instructions on the screen.

There are two download options: either download all logs or the last 10 only. During the loading of data in the USB memory, information about % of loading appears.

Press the **BACK** key to return to the settings menu or the **HOME** key to return to the initial screen.

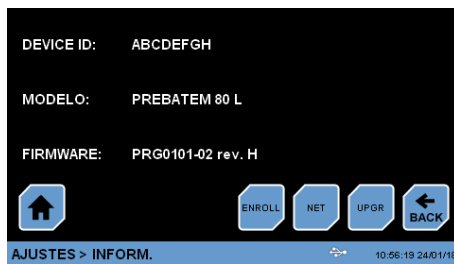
Fig. 8.9 Information screen

Fig. 8.9.1 Firmware update

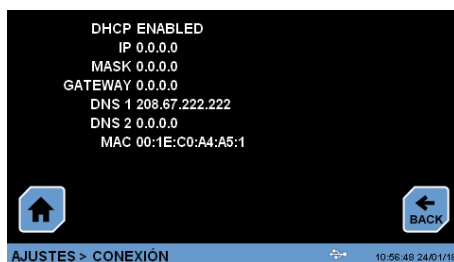
Fig. 8.10 Information about connection to the data network

8.9 Information. Firmware update.

The equipment is tested and checked before shipment but there may be improvements in the control operation.

If it is necessary to update the firmware version, press the **INFORMATION** option and Fig 8.9 appears.

Press the **UPGR** key and follow the instructions that appear on the screen. A screen similar to Fig 8.9.1 will appear.

To obtain the file with the update, it is necessary to know the **DEVICE ID** code that appears in screen Fig 8.9. Load this file into a **USB** memory and insert it into the front connector.

Attention: during the update process **DO NOT** turn off the equipment or remove the **USB** memory from the connector. The equipment could run out of control.

The update process lasts a few minutes. During the update process the control is turned off several times, but this is normal.

Press the **BACK** key to return to the settings menu or the **HOME** key to return to the initial screen.

8.10 Information about connection to data network

From the **INFORMATION** menu you can query communication data to a data network.

Press the **NET** button to see this information.

Press the **BACK** key to return to the settings menu or the **HOME** key to return to the initial screen.

Notice to customers:

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.