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# \*English Manual

## 1. Introduction

The S500 series of GSM temperature data logger is the new generation of data loggers; which has been added with the GSM alarm message function based on the original S500 series. Users can dial the data logger through any mobile phone to get the temperature value, besides, the temperature data logger will send alarm message to defined users automatically once the setting limits being exceeded.

### 1.1 Features

- Elegant appearance, easy to operate and reliable. Desktop and wall-mounted.
- High accuracy (0.2~0.5°C/2~5%RH) with large capacity. (65000 logs).
- Large LCD display, where temperature is shown.
- Internal audible and visual alarm and send alarm message to mobile phone once temperature is over limit.
- Data query function via mobile phone. Get the real time data anytime at anywhere.

### 1.2 Technical Specifications

Part number	2101520 / 2101522 (double probe)
Accuracy	±0.5°C
Measuring range	-40 to 85°C
Sensor type	Single temperature external sensor with 3 meters / 2101522 with double temperature probe
Capacity	65,000

### 1.3 Applications

It has been widely applied in agricultural research industry, food industry, medical industry, electrical industry, environment protection and laboratory fields.

### 1.4 Installation and instructions for SIM card

1. We suggest buying GSM SIM card (SIM card without monthly rent). It must be GSM network SIM card, not a CDMA network which is invalid.
2. Install or remove SIM card should be under the state of OFF in case any SIM identification failure and damage.
3. Just call the alarm temperature data logger via any mobile phone and it will respond you with a message when you want to know the temperature value.
4. Before installing the SIM card, you must ensure that you do not request the PIN code when starting it. The computer is unable to communicate.


## 1.5 Notice

- 1、Only GSM network SIM card is available.
- 2、Ensure that the SIM card is of certain amount of money for the text message fee.
- 3、The data logger must be supported by External Adapter—DC 9V and turn on the alarm function to activate the alarm message system, otherwise it cannot send the message successfully.**

## 1.6 Instructions


1. Be sure that the battery is installed, external power connected and SIM card insert well before you turn on the data logger.
2. Install the software and connect the data logger with PC via USB connection cable.
3. Run the software and click the “connect” button. The factory communication settings between the PC software and the data logger are already used to communicate the two devices. If necessary, modify the parameters of the PC software to adjust.
4. Click **【setup】** and set the high and low limits as well as the log interval. The name of the data logger should be in English and no illegal symbol. (Notice: Data logger will send alarm message when setting limits are exceeded and the data logger record the data. Suggest set log interval above 60 seconds.)
5. Click **【Options>GSM】** in the tool bar and set the mobile phone number. It can set 3 numbers in total.

## 1.7 Product structure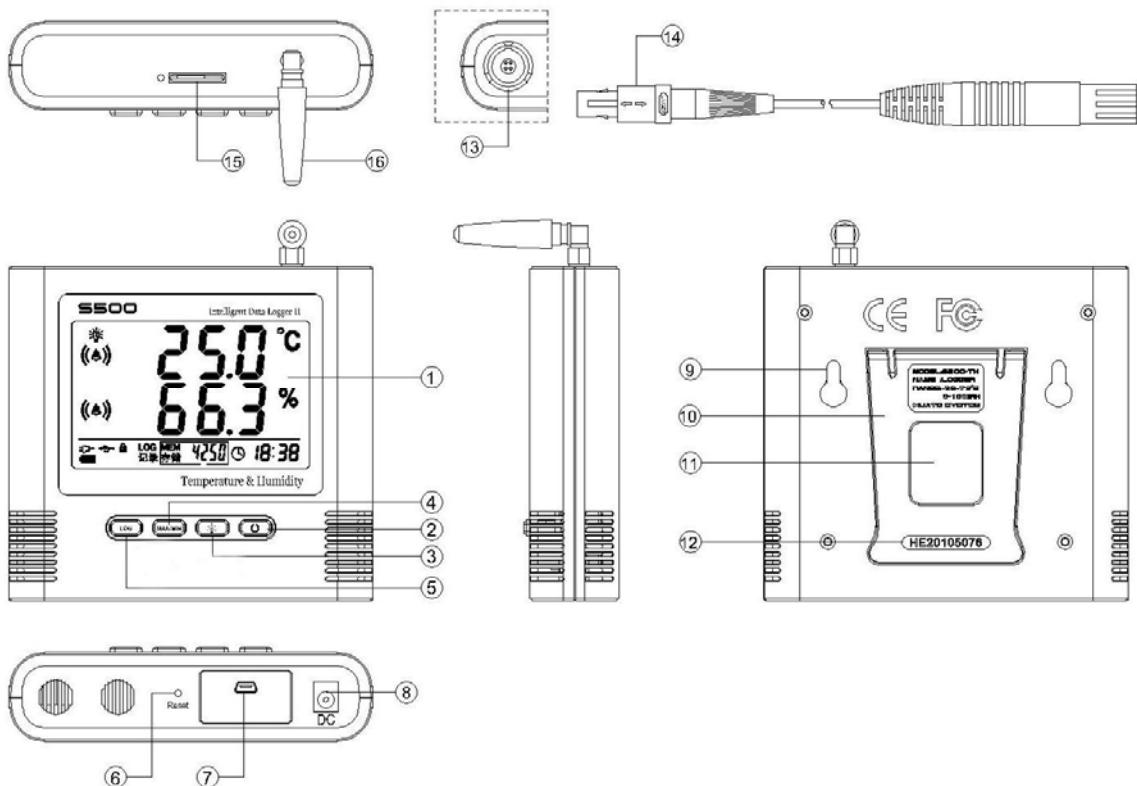

LCD display	9. Hanger (used for fixing on the wall)
ON/OFF	10. Holder (used for desktop)
Backlight	11. Battery cover (Screw to replace the battery)
MAX/MIN checking button	12. Serial number
LOG	13. Probe connector (two connectors on model 2101522)
Reset button	14. Temperature probe
USB port	15. SIM card slot
DC 9V port	16. Wireless antenna

## 1.8 LCD display instructions

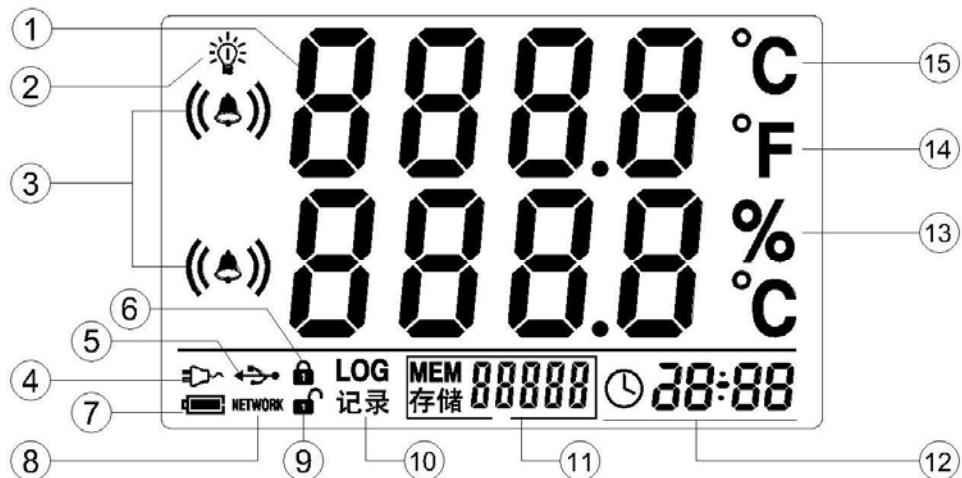- ① Temperature values display (double display on model 2101522)
- ② Symbol for backlight.
- ③ Symbol for alarm, it will alarm when temperature setting limits will be exceeded.
- ④ This symbol shows that data logger is being supported by external power—DC 9V.
- ⑤ This symbol shows that data logger is connected with PC and you can download all the recording data.
- ⑥ This symbol shows that all buttons are locked.
- ⑦ Battery indication
- ⑧ This symbol shows that SIM card inset well.
- ⑨ This symbol shows that buttons are unlocked
- ⑩ This symbol shows that data logger is being the recording status.
- ⑪ Display area for recording numbers
- ⑫ Display area for year/month/date/time; alternate display.
- ⑬ Symbol for humidity unit
- ⑭ Symbol for temperature unit
- ⑮ Symbol for temperature unit

## 1.9 Instructions for buttons

 LOG

Log function button, turn on/off the log function.

 MAX/MIN

Maximum and minimum button. Press to check the maximum and minimum value of all the recording data.

Backlight on/off button.

On/Off button

### Notice:

- 1) Press  and  simultaneously to lock and unlock data logger.
- 2) Data logger has alarm message system only when data logger is under logging status.

## 2. Software

### 2.1 Requirements of the Computer Hardware

1. OS: Windows XP/Vista/7/8 (32/64bit)
2. CPU: 1.6GHz
3. Physical Memory: 512MB
4. Hard-drive Space: 4 GB
5. USB Port: 1

### 2.2 Install Driver & Software

#### 1. Install Driver:

Notice: If it is Windows 64bit, please install 64bit-driver.

#### 2. Install Logpro Software:


To install the software, simply copy the files to a known folder. We recommend creating a link to the desktop to have direct access to the .EXE file.

### 2.3 Setting the logger's attribute to start a new measurement.


1. Connect the data logger to the computer.<sup>1</sup>
2. Running Logpro software on the PC.
3. From the toolbar select Connect.<sup>2</sup>

<sup>1</sup> Windows operating system cannot handle USB devices being unplugged and plugged back too fast. When unplugging the logger, wait for about 5 seconds before plugging it in again. If you unplug and plug back a device too quickly, the computer may stop recognizing any USB devices on that port. If this happens, you will have to restart the computer. This is a windows USB problem and is not related to Logpro.



<sup>2</sup> If more than one logger is connected to your computer at the same time, the program will ask you to choose COM port manually.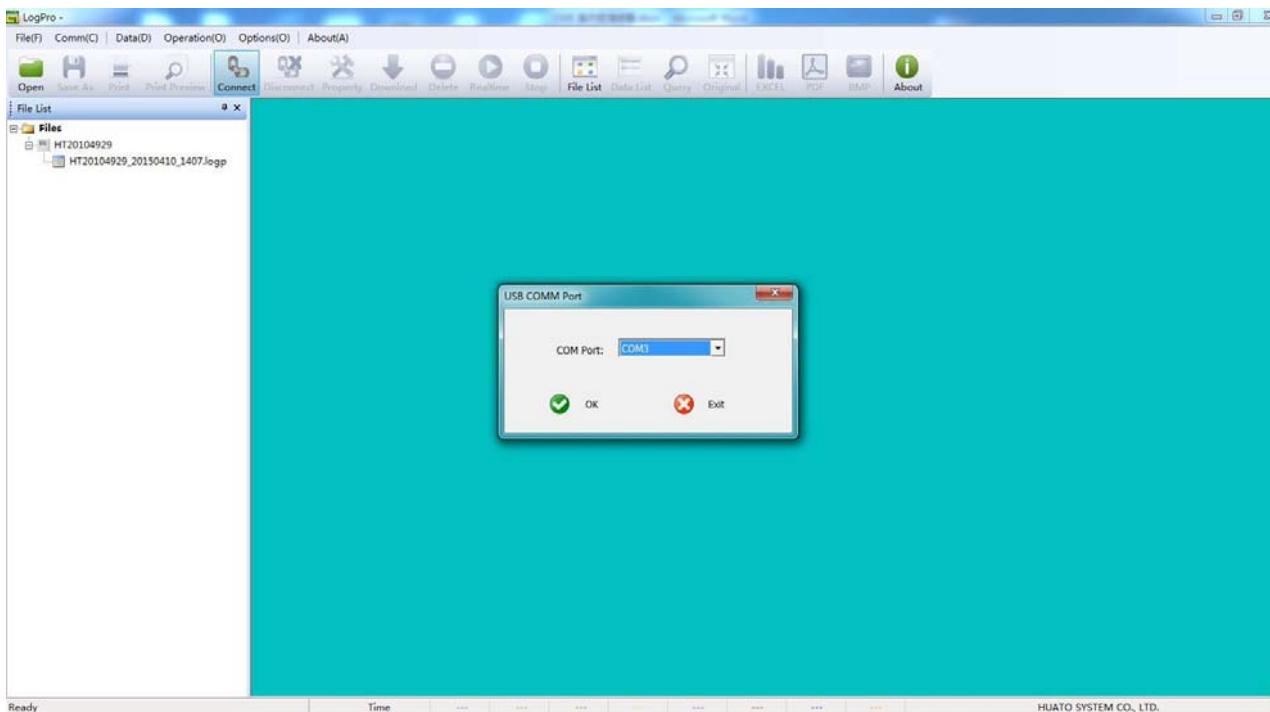


4. Read the logger's attribute.
5. Set the attribute.
6. Sync the attribute & time. (The PC time will be synchronized to the logger as well.)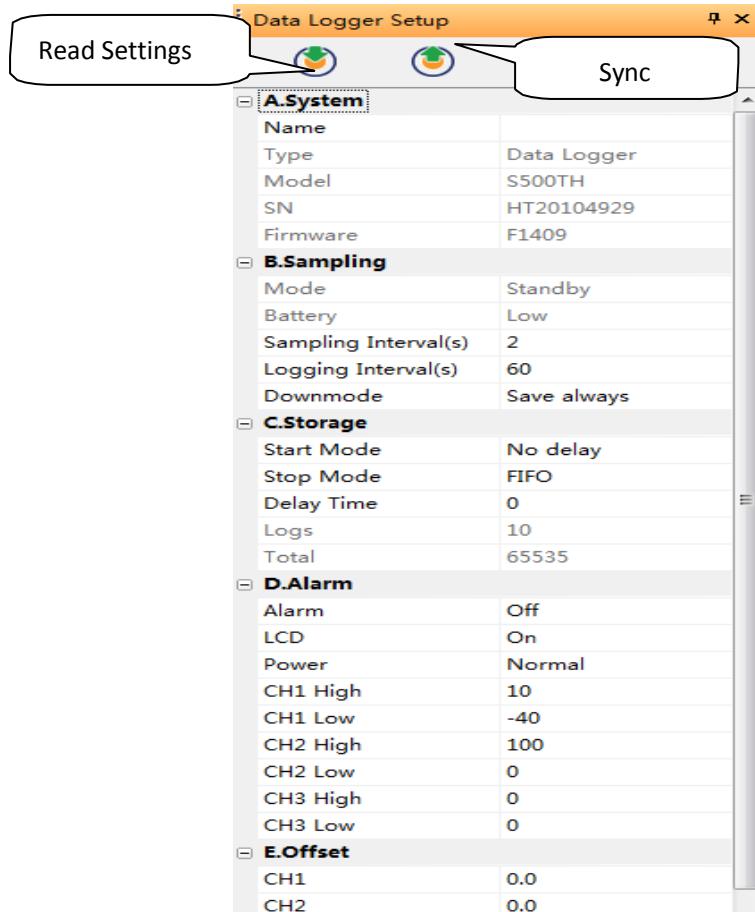

7. Choose (**GSM setup**) to setup cell phone number to receive alarming message.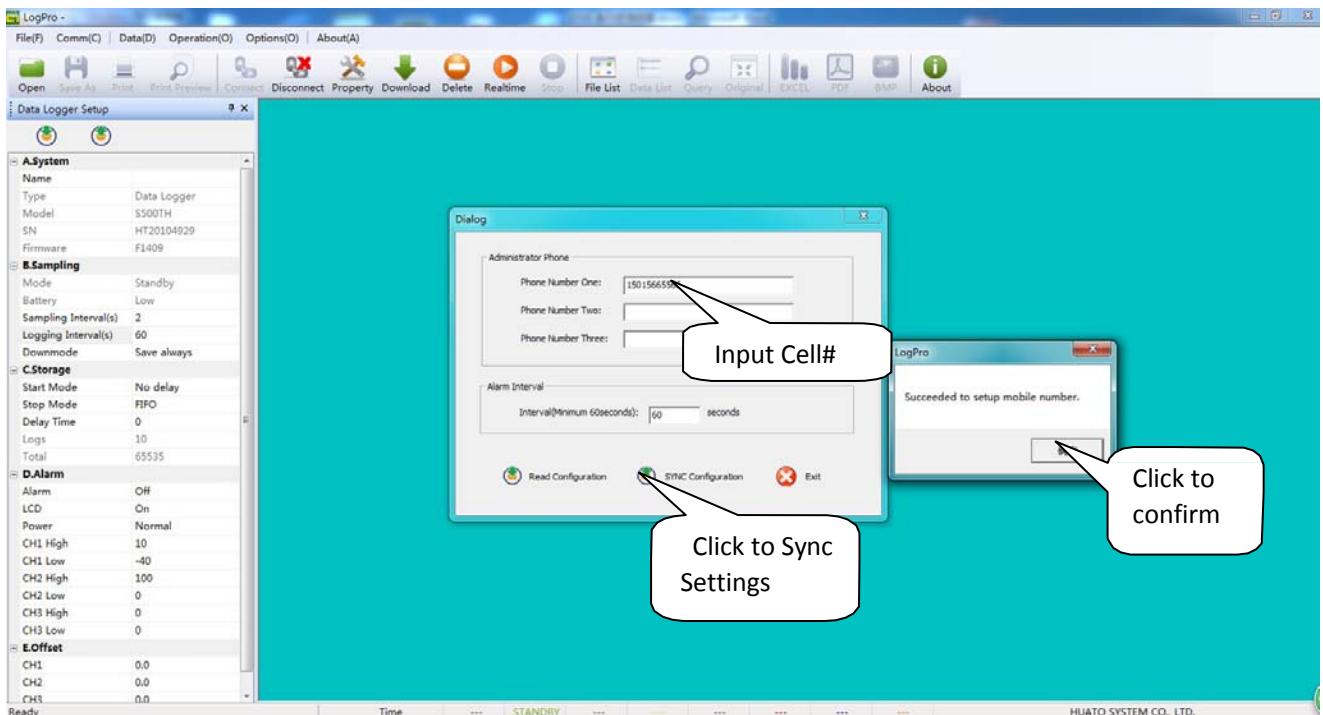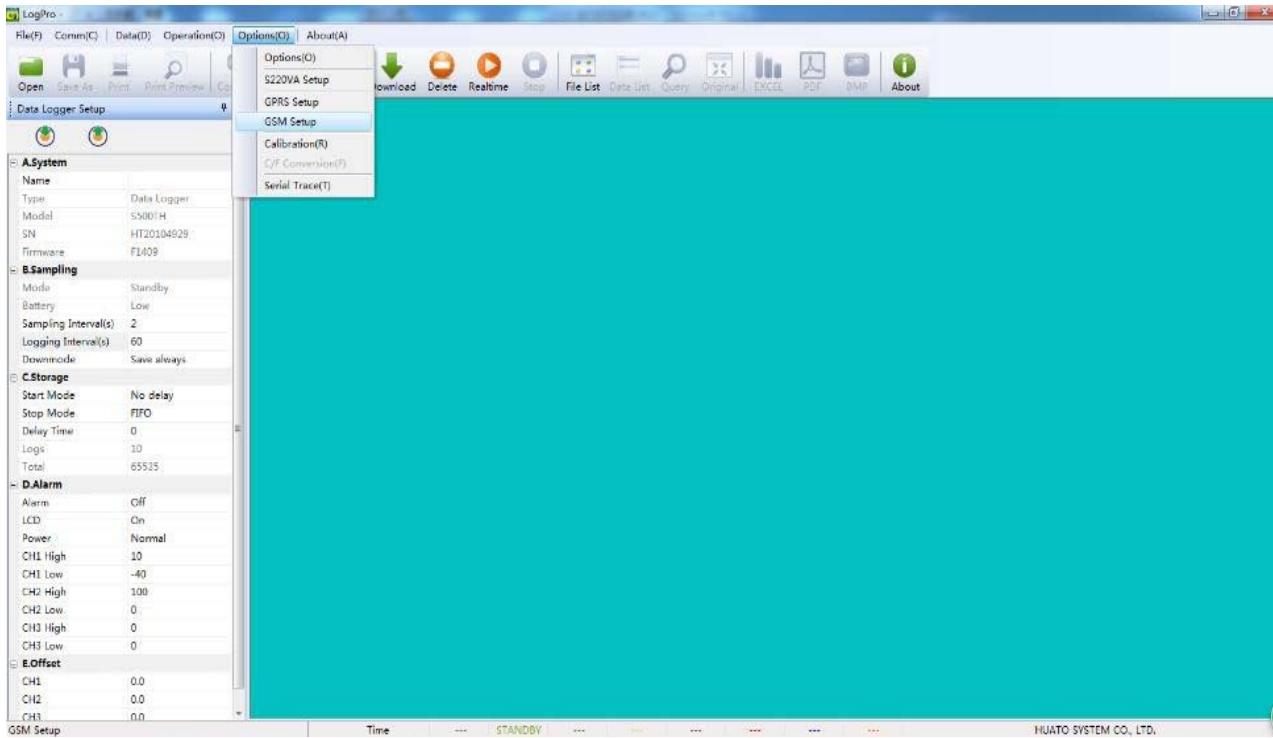

## 8. Setup the temperature and humidity upper and lower limits.

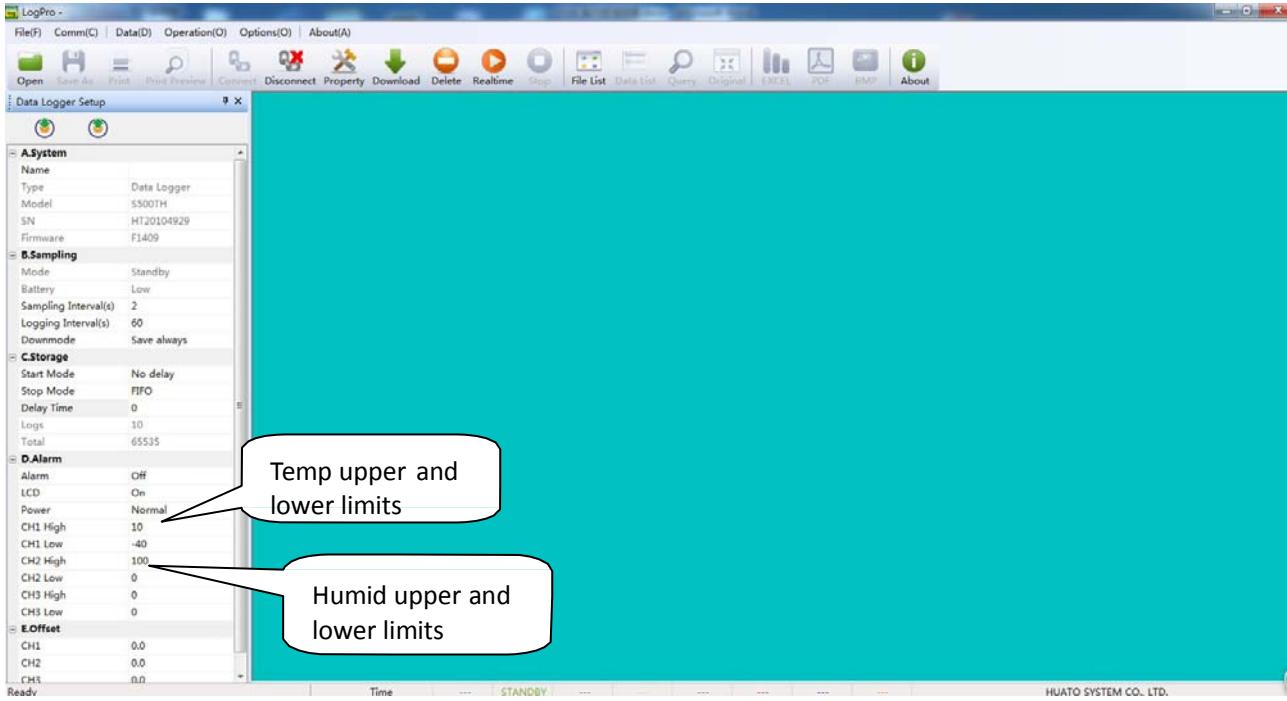
## 9. Then sync the setup with loggers after finish setup.


## 2.4 Data Analysis

Before starting, set the folder where to save the data inside **Options> Options**.

1. Click on (**Download**) to download data to the computer, it will be automatically saved and will generate reports and data list.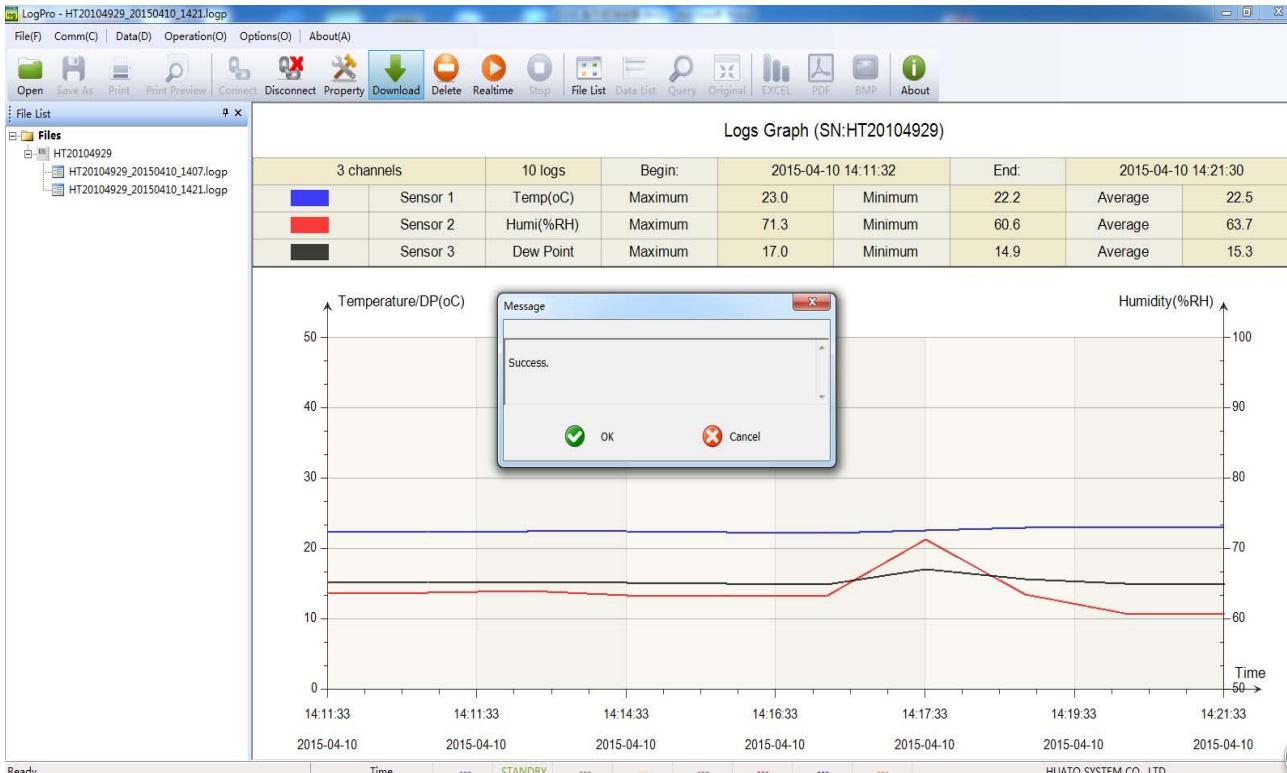

2. Click on (**Data List**) to display downloaded data on the left.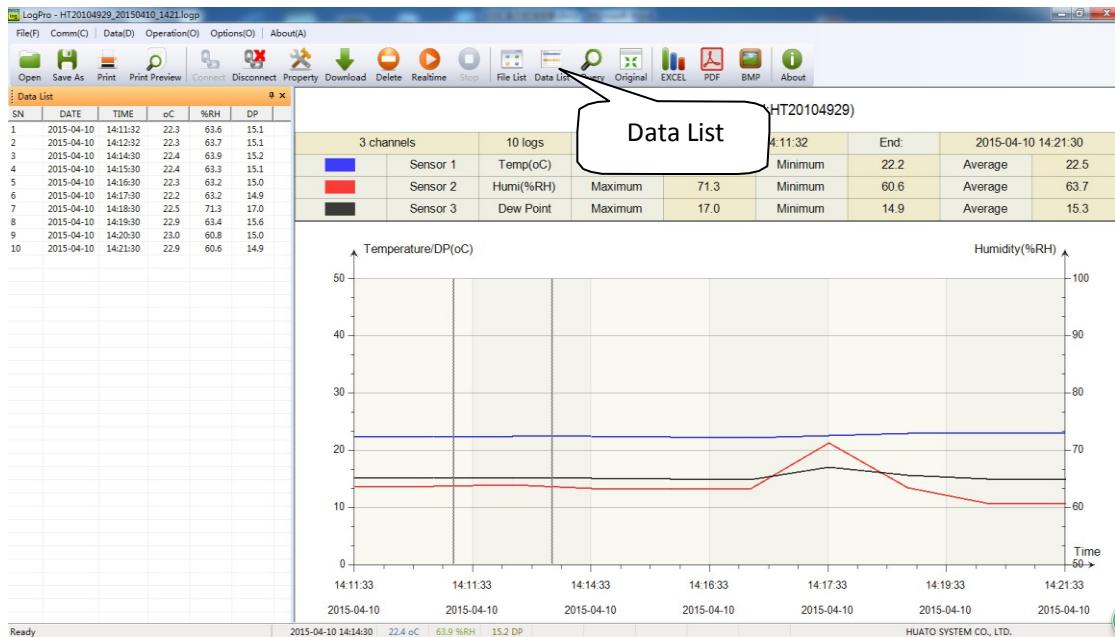

3. Click on to save the data in the Excel format; or click on to save the data in a PDF

file; or click on to save data in a picture.

4. Click on (**Print**) to print data or (**Delete**) to delete data.