**REACTOR PARA PLANTA PILOTO
REACTOR FOR PILOT PLANT**

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INDEX

1. General	35
1.1 Introduction	35
1.2 Safety.....	35
1.3 Instrument Synopsis	38
1.4 Instrument Configuration.....	39
2. Unpacking and Installation	40
2.1 Prepare	40
2.2 Installation.....	40
2.3 Reactor and associated devices connection.....	47
2.4 Operation.....	48
3. Technical Data	49
3.1 Borosilicate Glassware 3.3	49
3.2 Lubrication.....	50
<i>Notice to customers</i>	50
4. Operating	51
4.1 Caution	51
4.2 Operating.....	52
5. Maintenance and Management	53
5.1 Management	53
5.2 Cleaning.....	53
5.3 Maintenance	53
5.4 Troubleshooting	54
6. Accesories and Spare Parts	59
7. Examples of Use of Support Trolleys	64
8. Examples of Application	65

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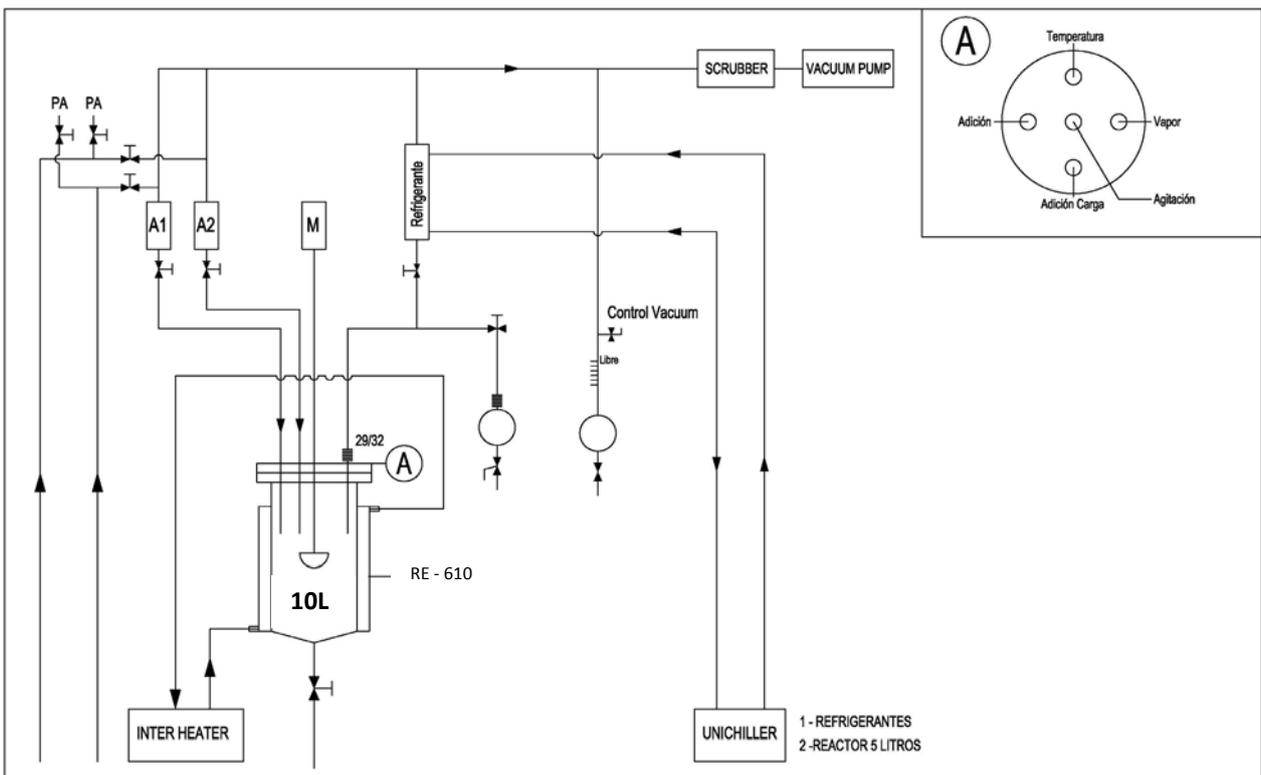
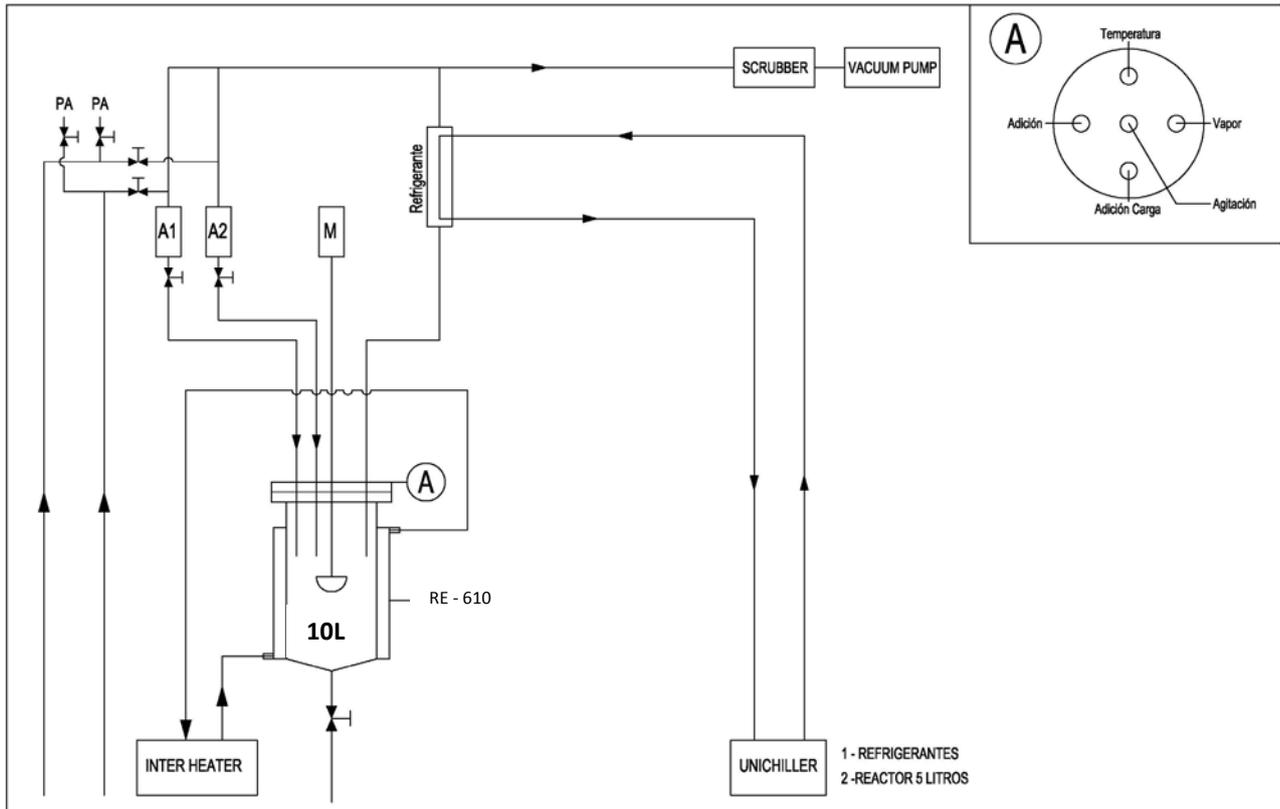
1. General

1.1 Introduction

Thanks for using 'GR-series speed-regulation glass reactor'. This operation manual stated the installation, using, maintaining and servicing of glass reactor. Before using this instrument, please ensure that you have read and understood this manual completely.

1.2 Safety

This chapter describes the installation, safety rules in the process of using of 'GR-series speed-regulation glass reactor'. Users must grasp the related warning signs, strictly abide by the operation procedures to ensure the security of the equipment and personal and avoid the occurrence of accident.

1.2.1 User's Qualification

GR-series speed-regulation glass reactor must be operated by the person who has the practical operating experiences and can grasp of the detailed requirements in this manual. Otherwise, it must be used under the guidance of the person who has the related technology skills.

1.2.2 Proper Use

Application fields:

- a) The process which need high temperature, low temperature conditions.
- b) The process which need reduce pressure and stirring at the same time.
- c) Dispersion process.
- d) Distillation (reflux), concentration.
- e) Assembled into a glass reactor total system according to the specific requirements.

1.2.3 Improper Use

The operation that not abide by the related stipulation in this manual are regarded as improper use. Any damage caused by improper use is responsible by the users themselves.

Operation under the following conditions is prohibited:

- a) Used in explosive gas environment or explosive dust environment.
- b) Used in the places which the power supply is not in conformity with the requirements.
- c) Used in high magnetic fields, corrosive environment.
- d) Hard, brittle bulk materials (such as stones, soil samples, etc.) may damage the vessel.
- e) The volume of the sample added to the vessel exceeds the specified limit.

1.2.4 Warning Sign

Sign	Description
 <p>Danger</p>	<p>Danger It shows that the situation is very dangerous and will lead to death or serious injury.</p>
 <p>Warning</p>	<p>Warning It shows that the situation is very dangerous and will lead to death or serious injury.</p>
 <p>Prohibit</p>	<p>Prohibit Get wet in the rain or splashing water.</p>
 <p>Note</p>	<p>Note It shows that the situation is very dangerous and will lead to injury.</p>
	<p>Note Caution high temperature.</p>
	<p>Note Please wear protective equipment, otherwise may cause personal injury.</p>
<p>Reminder</p>	<p>Reminder May cause equipment damage.</p>

Especially pay attention to each warning sign.

1.2.5 Hazards Related to the Instrument

Please pay attention to the following safety tips:

 <p>Danger</p>	<p>Danger</p> <ul style="list-style-type: none"> ● Do not use it in explosive gas environment or explosive dust environment.
 <p>Warning</p>	<p>Warning</p> <ul style="list-style-type: none"> ● Make sure the power supply is in conformity with the requirements on the name plate ● Equipotential connection has been done, do not remove inner grounding connection line. ● Before open the equipment enclosure for maintenance and repair, be sure to disconnect the power supply, then operate it after 5s, ensure the residual voltage release to safety value to avoid electric shock. ● Do not use broken glass parts
 <p>Prohibit</p>	<p>Prohibited</p> <ul style="list-style-type: none"> ● Do not use it outdoor. Get wet in the rain or splashing water will cause electrification on metallic shell surfaces, and will cause casualties.
	<p>Warning</p> <ul style="list-style-type: none"> ● Do not touch the vessel surface and connection pipes with your body directly to avoid scalding.
 <p>Note</p>	<p>Note</p> <ul style="list-style-type: none"> ● If there's any abnormal situation when using it, please disconnect the power supply immediately for troubleshooting or contact professional maintenance personnel. ● Be aware of broken glass parts. ● It may have static electricity when adding material to the vessel through feeding port
<p>Reminder</p>	<p>Reminder</p> <ul style="list-style-type: none"> ● Do not use it in the environment mentioned above. Otherwise, it will affect its normal running and lifetime. ● Keep well ventilated around the equipment. ● Do not use it when there's no one there.

1.2.6 Other Hazards

 <p>Warning</p>	<ul style="list-style-type: none"> ● The sample in the vessel or some solvents near it may form peroxides, high concentrations of flammable solvents, etc. ● Acid, alkali steam and other corrosive materials around the equipment will damage the equipment insulation and affecting component performance and service life. ● Beware of the risk of explosion when dealing with hazardous materials or unknown samples.
--	--

1.2.7 Safety Measures

	<ul style="list-style-type: none"> ● Please wear personal protective equipment when operating this instrument, such as protective glasses, protective clothing and gloves
---	--

1.3 Instrument Synopsis

RE-610 speed-regulation glass reactors are mainly used for synthetic reaction, distillation and concentration of different types of materials. We have single wall, double wall and triple wall glass reactors.

The reactor can be pumped to a negative pressure state according to some experimental requirements. A constant pressure funnel or an adjusting valve on the feeding bottle allows you to add material the vessel at a uniform and controlled speed. Condenser helps to recover some materials distilled during reaction.

Single wall glass reactor has no jacket.

Double wall glass reactor has a jacket. Circulating liquid can be filled into the jacket to heat or cool materials inside.

Triple wall glass reactor has two jackets, the inner jacket is for circulating liquid, and the outer layer should be pumped to vacuum condition for thermal insulation.

1.4 Instrument Configuration

RE-610 speed-regulation glass reactor (Customized product as per the real object)

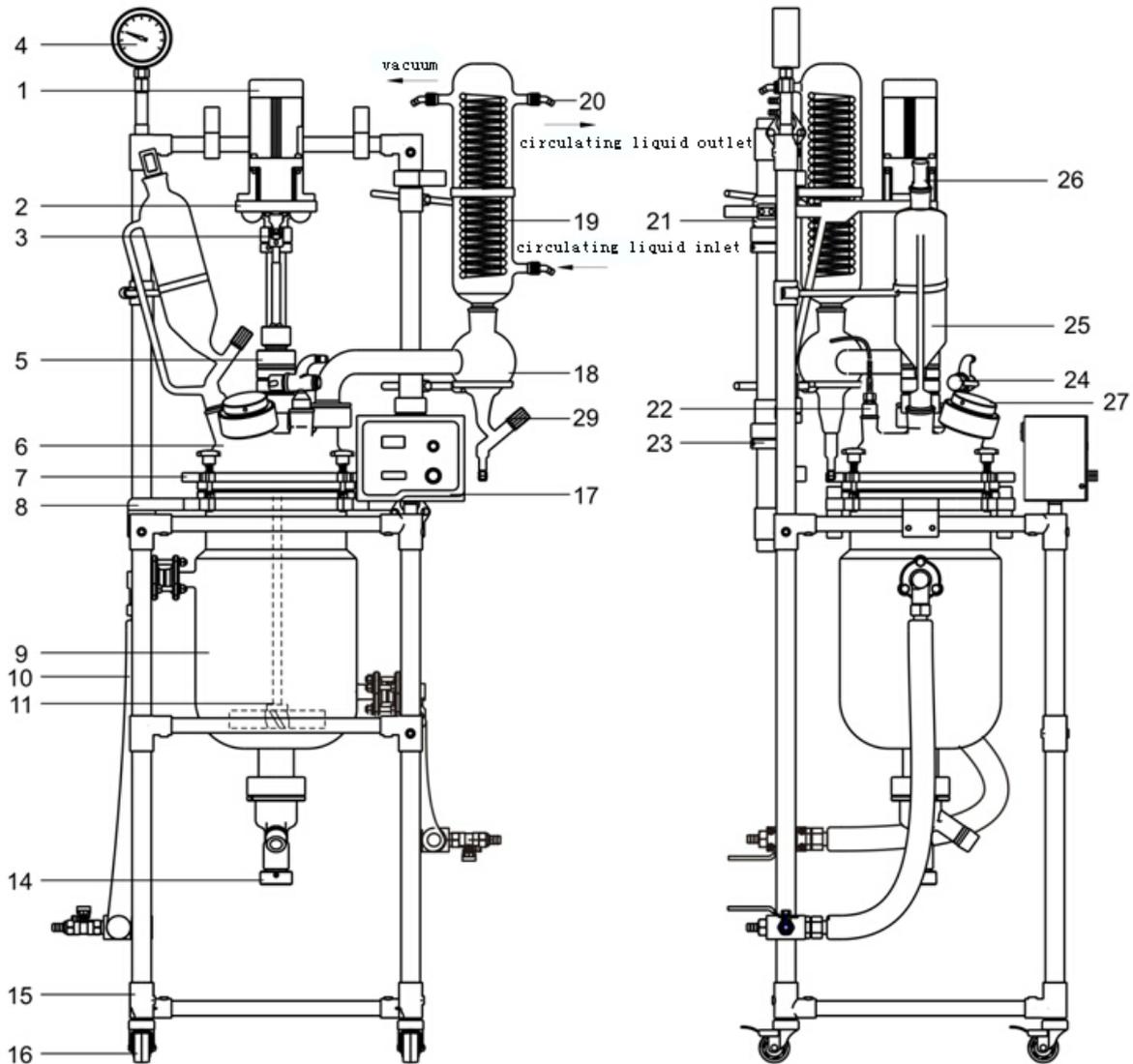
Figure 1.4.2 - RE-610 Speed-regulation glass reactor

1 Motor	2 Motor tray	3 Universal coupling	4 Vacuum gauge
5 Stirrer bearing	6 Reactor lid	7 Upper tray	8 Middle tray
9 Vessel	10 Circulating pipe kit	11 Stirring paddle	14 Discharge valve
15 Frame	16 Caster	17 Control box	18 Condensing dispenser
19 Condenser	20 Joint	21 Fixing frame	22 Sensor
23 Fine tuning unit	24 Feeding valve	25 Constant-pressure funnel	26 Plug
27 Solid charging port sealing kit	29 15# cock		

2. Unpacking and Installation

2.1 Prepare

1) After unpacking the case, please read the operation manual carefully, and check whether the components are enough compare with the packing list.

2) Take out the frame, put it on the floor then lock the casters.

Note: The vessel and frame are assembled completely of reactors less than 20L.

3) Please clean the glass parts to maintain its cleanliness to meet the test requirements before installation.

Note: Glass device is fragile, please be careful when unpacking.

2.2 Installation

2.2.1 RE-610 vessel installation

1) The vessel (9) fits from the bottom of the middle tray (8).

2) The sealing ring (71) is fitted from the flange mouth of the vessel (9) .

3) Put the sealing ring (71) on the middle tray (8) to make their circle fitted.

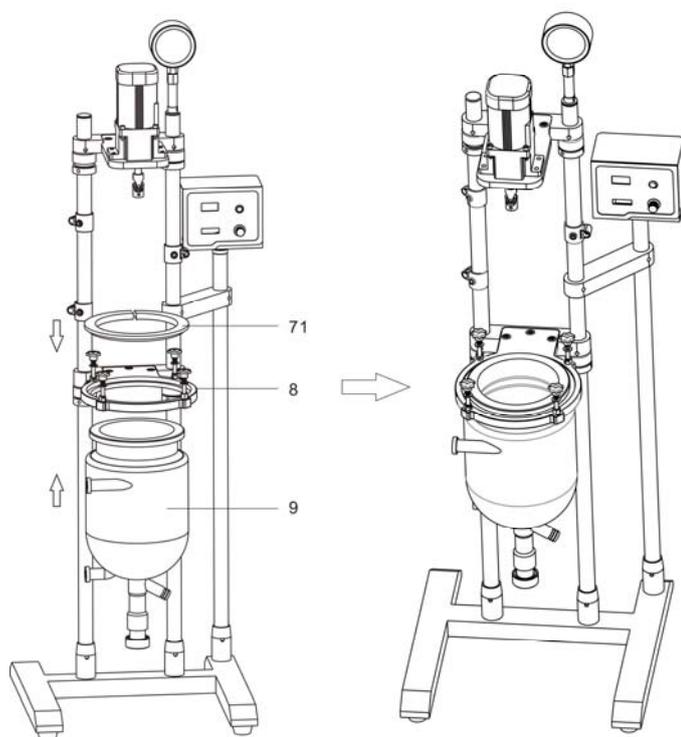
Figure 2.2.1 Vessel installation of RE-610

2.2.2 Stirring paddle installation

Four plum knobs loose to the bolts can be flat, remove the upper tray (7), tray gasket, reactor lid (6) and reactor lid openings sealing gasket. Place the stirring paddle (11) in the vessel body. Coated vacuum grease on the contact surface of the vessel body and the reactor lid. Install the reactor openings sealing gasket, reactor lid (6), tray gasket (7) and upper tray (7) in turn, tighten the plum knob.

Note: Please refer to 1.4 for the mounting position of the reactor lid.

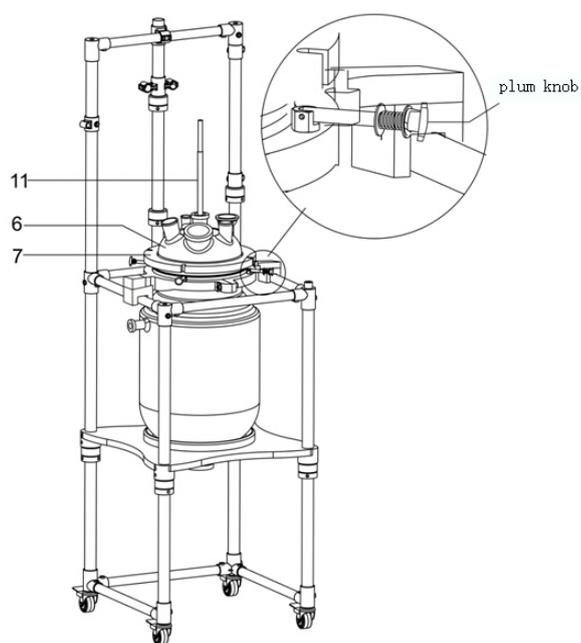
Figure 2.2.2 Stirring paddle installation

6 Reactor lid 7 Upper tray 11 Stirring paddle

2.2.3 Stirring shaft installation

Please install the stirring paddle according to figure 2.2.3.

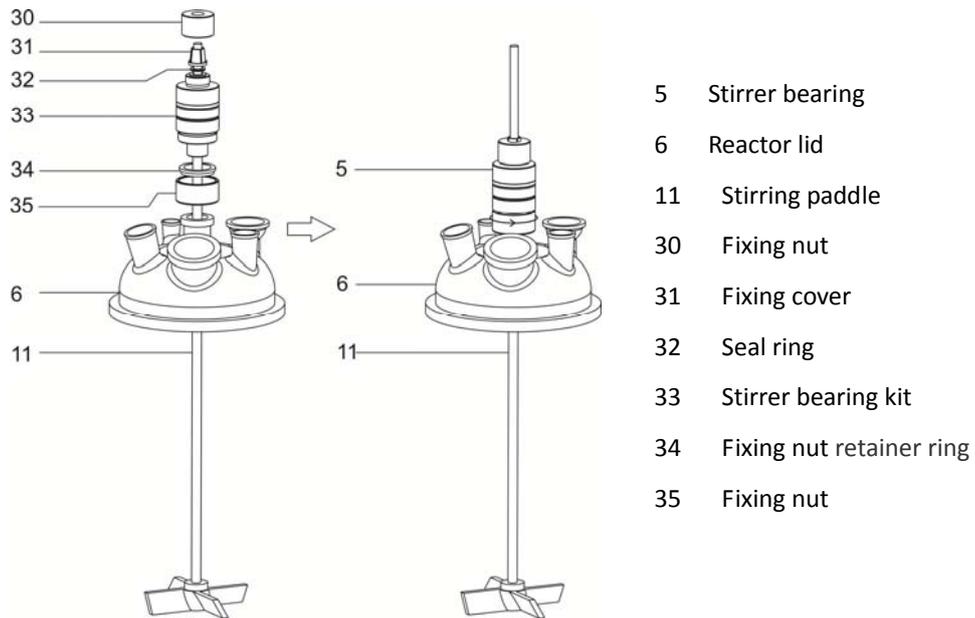
Figure 2.2.3 Stirrer bearing installation

2.2.4 Motor installation

The motor, motor tray, universal coupling and fixing pole have been installed together before leaving factory, please see figure 2.2.4.1.

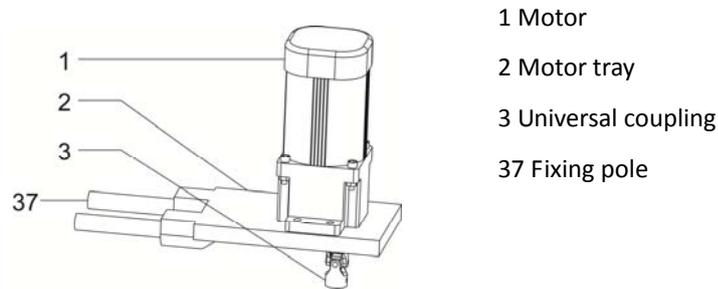
Figure 2.2.4.1 Motor assembly

1) RE-610 motor installation

Install the motor kit on the fixing frame (21) to make the motor axis and stirring paddle coaxial, tighten universal coupling (3). Adjust the height of the motor fixing frame (21) to make the gap between the lower of the stirring paddle and the vessel bottom >30mm, fix the fixing frame (21). Tighten the fixing nut (30) to fix the stirring paddle.

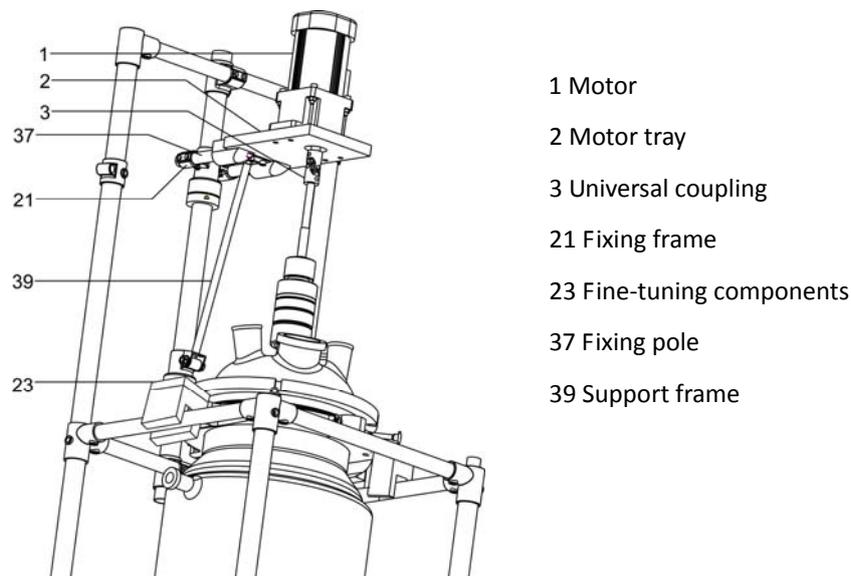
Figure 2.2.4.3 RE-610 motor installation

Connect support frame (39) and motor tray (2), tighten the connection screw. Install the motor kit on the fixing frame (21) to make the motor axis and stirring paddle coaxial, fix the universal coupling (3). Adjust the position of fixing frame (21) to make the gap between bottom of the stirring paddle and vessel bottom >30mm, fix the fixing frame (21). Then tighten the connecting bolt and screw on the lower position of the support frame (39). Adjust the fine-tuning components (23) to make the level of the motor tray is horizontal. Tighten the fixing nut (30), fix the stirring paddle.

2.2.5 Nozzle and condenser installation

Please install it according to figure 2.2.5.

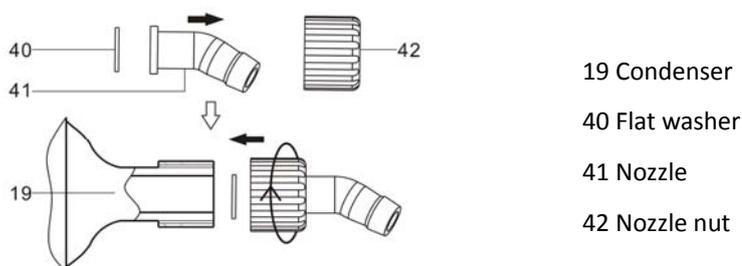
Figure 2.2.5 Nozzle and condenser installation

2.2.6 Condensing dispenser and condenser installation

RE-610 Condensing dispenser and condenser installation

Place the tray (44) on the frame, put the condensing dispenser (18) on the tray (44). Connect the condensing dispenser (18) and the reactor lid joint, adjust the position of tray (44), tighten the fixing screw (45) and fixing nut (48). Put the condenser (19) on the upper joint of the condensing dispenser (18), the rubber bracket (43) is locked around the upper portion of the condenser (19), please see figure 2.2.6.2.

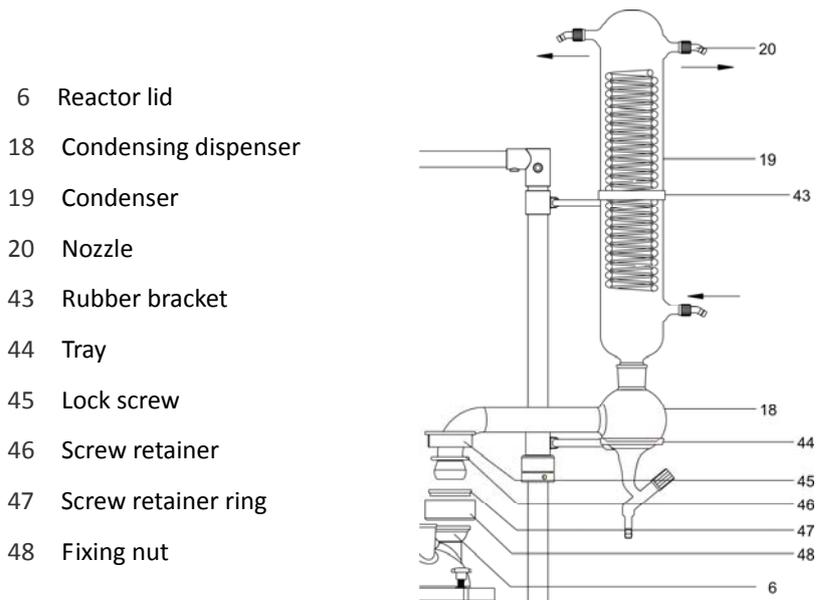
Figure 2.2.6.2 RE-610 Condensing unit installation

2.2.7 Sensor installation

Put the fixing screw (49), fixing cover (50), O-ring (51) and fixing base (70) on the sensor in turn, adjust sensor (22) position, tighten fixing nut (49).

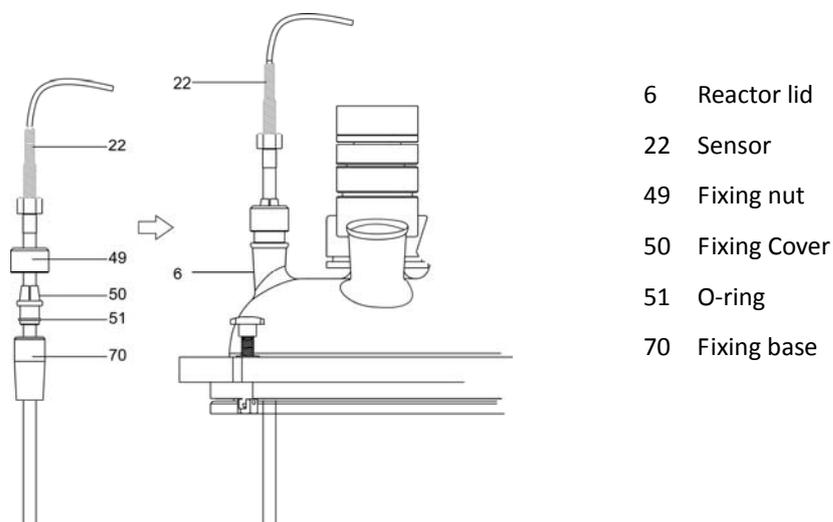
Figure 2.2.7 Sensor installation

2.2.8 Solid charging port sealing kit installation

Please install it according to figure 2.2.8.

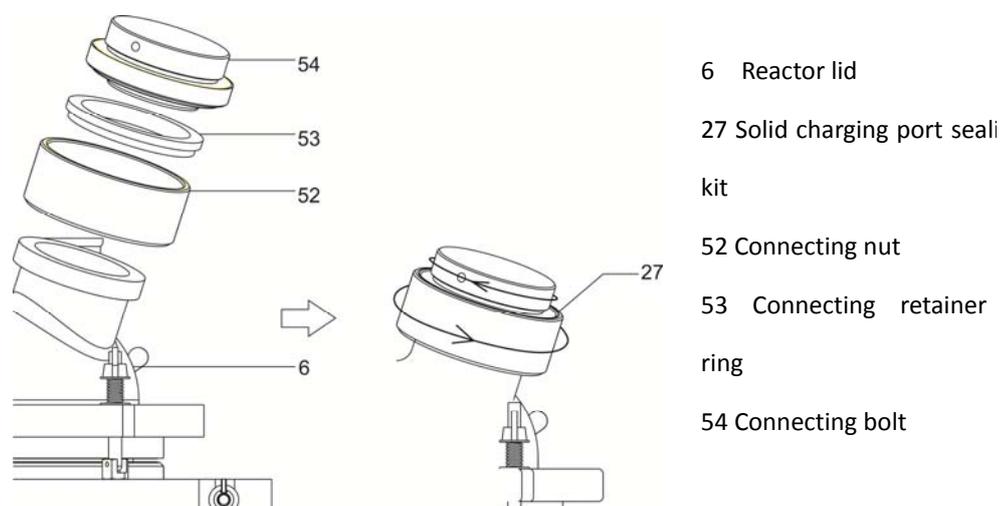
Figure 2.2.8 Solid charging port sealing kit installation

2.2.9 Constant-pressure funnel and feeding port valve installation

Please install it according to figure 2.2.9.

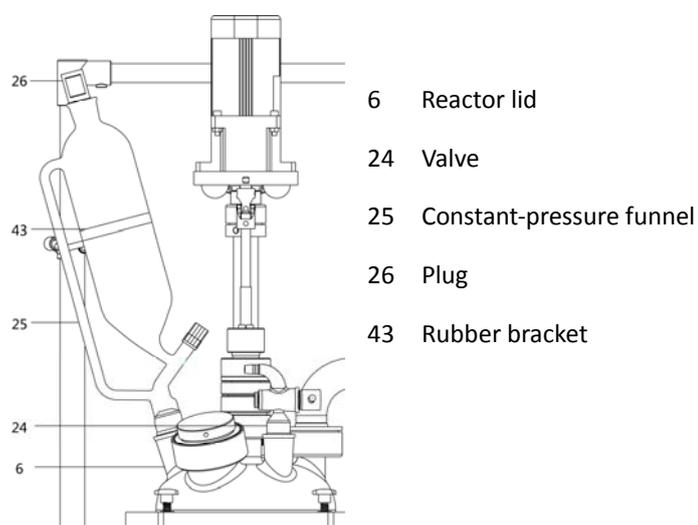
Figure 2.2.9 Constant-pressure funnel and valve installation

2.2.10 Discharge valve installation

Note: Please open the valve (14) before install it in case of damage the vessel body.

1) RE-610 discharge valve (14) installation

Put 80# flange mouth screw (59) and 80# flange mouth retainer ring (60) on the flange mouth of vessel body (9) and discharge valve (14), put 80# flange mouth sealing ring (61) on the flange mouth of discharge valve (14), position it in place, tighten the butterfly nut.

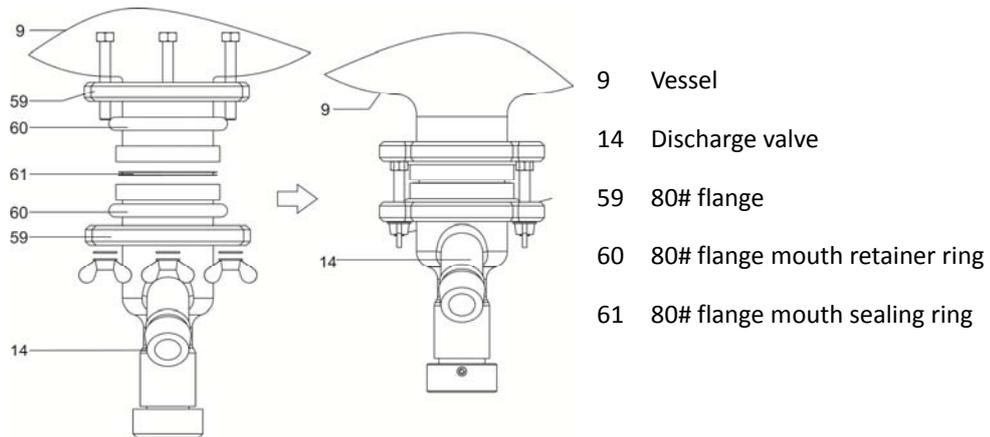
Figure 2.2.10.1 RE-610 discharge valve installation

2.2.11 Vacuum gauge and control box installation

Please see chapter 1.4 to find the installation position of Vacuum gauge (4) and control box (17).

2.2.12 Vessel jacket circulating pipe kit installation

The lower connecting joint of the vessel jacket is inlet, the upper is outlet. Please refer to figure 2.2.10.2 to install it.

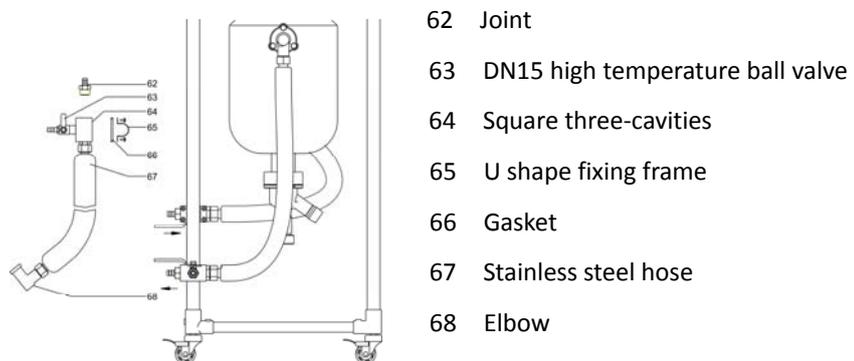
Figure 2.2.12 Circulating pipe kit installation

2.3 Reactor and associated devices connection

Connect the reactor with Dynamic Temperature Control System and vacuum pump, make sure that there's no leakage between each joint.

Note: One of a nozzle of the vacuum gauge connected with the upper vacuum joint on the condenser, and another nozzle connected with vacuum pump.

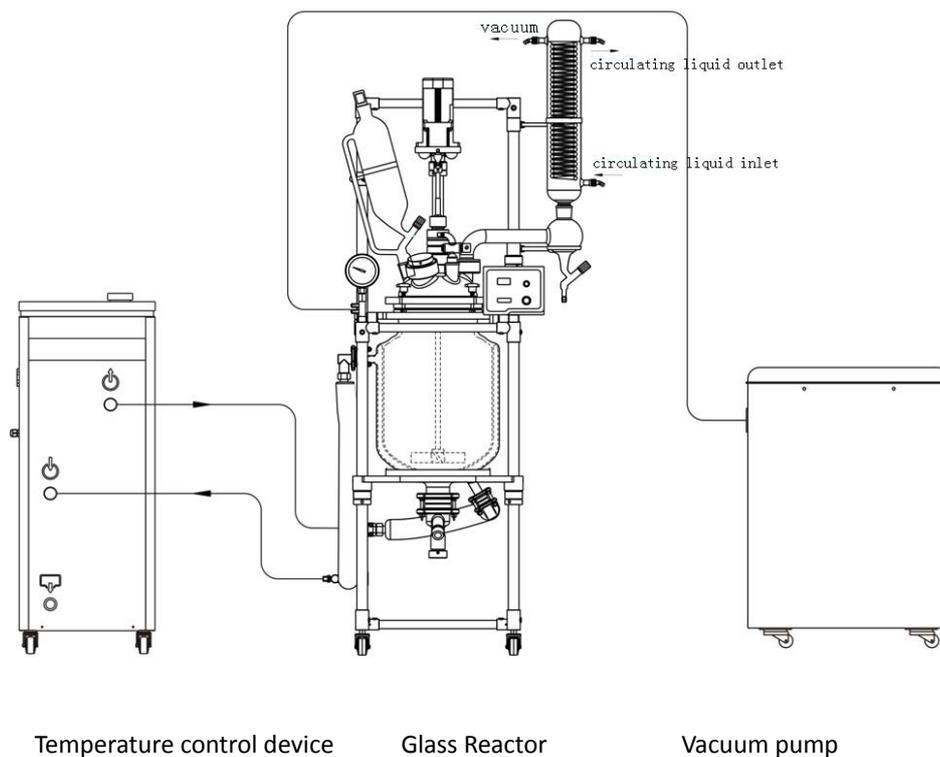
Figure 2.3 Total system connection

2.4 Operation

1. Turn the power switch on and the temperature and the speed will be displayed.
2. The speed is adjusted to 0, click on the Rotation key, slowly adjust the speed, the rotation of the agitation blade, and observe if any anomaly occurs.
3. Pour the materials into the container through solid feed or other inputs.
4. Evacuate the container according to the reaction needs. For the triple reaction vessel, the vacuum pump can be applied to evacuate the outer layer, to achieve the thermal insulation functionality.
5. Set the temperature control device according to the temperature shown, to obtain the desired reaction temperature inside the container.
6. Slowly adjust the knob to achieve the desired rotation speed.
7. Let the water or coolant pass through the condenser according to the process requirements.
8. After the reaction, first turn the heating or cooling temperature control device function off, maintain the cycle and then close the circulation function when it is near to the ambient temperature.

***Note:** The temperature control unit cycle is turned off. The circulation valve can not be closed, leave it open. Otherwise, if the body of the container is too hot or too cold, it could cause inflation and cause damages to the equipment.

9. After the speed test is set to 0, click on the key 1 minute after closing the reactor power
10. Open the discharge valve to discharge the materials. If necessary, the valve parts or the sealing parts of the container should be cleaned and then reassembled safely.

***Note:** The glass parts must be cleaned after returning to normal temperature. The droplets on the inside and outside should be cleaned after washing to avoid rupture of the container, caused by an irregular temperature or by a sudden change in temperature.

3. Technical Data

Table 3 - RE-610 speed-regulation glass reactor technical data

Model	RE-610
Stirring speed (rpm)	50~500
Jacket pressure (Mpa)	Inlet≤0.03, outlet 0
Pressure in the vessel (Mpa)	-0.1~0.0
Bearable temperature difference (°C)	-80~200
Motor rated power (W)	90
Glass vessel capacity (L)	10
Jacket capacity (L)	About 3
Vessel jacket heat exchanging area (m ²)	0.25
Condenser heat exchanging area (m ²)	0.16
Diameter of circulating fluid inlet and outlet	DN15
Reactor lid openings	6
Constant-pressure funnel (L)	1
Receiving flask (L)	--
Power supply	220 - 240V~, 50/60Hz
Operating temperature (°C)	5~35
Relative humidity (%)	≤ 65
Enclosure protection grade	IP20
Dimensions (mm) W×D×H	625×820×1780
Drain port ground clearance (mm)	About 355
Net weight / Gross weight (kg)	50/90

3.1 Borosilicate Glassware 3.3

- Chemical properties
- Hydrolytic resistance
- Acid resistance
- Alkali resistance
- Temperature resistance
- Optical

Technical features of working use in standard reactors: RE-610

Vacuum: -0.9 to +0.5 pressure bars

Working temperature: -30°C to +200°C

PTFE gaskets: -100°C a +250°C (maximum working -20 to +200°C)

Glass is an organic mixture of metal oxides fused together at high temperatures, which upon cooling, solidifies into the clear, rigid, non-crystalline, versatile material known widely across the globe

Borosilicate glass has a very high resistance to attack from water, acids, salt solutions, halogens and organic solvents. Only hydrofluoric acid, hot concentrated phosphoric acid and strong alkaline solutions cause appreciable corrosion of the glass.

Borosilicate glass has excellent thermal properties at both high and low temperatures. The maximum recommended working temperature for laboratory glassware manufactured is 500°C (for short periods of time only). Special care should be taken at temperatures above 150°C to ensure both heating and cooling is achieved in a slow and uniform manner.

3.2 Lubrication

High vacuum grease

Range Hg: 10-6 mm

Range: -40°C to 200°C invaluable

This grease does not melt and is been composed for lubrication and sealing of glass and wrenches in high vacuum systems with outstanding stability.

Notice to customers:

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.

4. Operating

4.1 Caution

Table 4.1 Caution

 <p>Note</p>	Please read this manual carefully before using it.
	Put it on the floor then lock the casters.
	The best ambient temperature for this device is 5~35°C. It will be broken if working outside this temperature range for a long period of time.
	Avoid strong vibration, otherwise it may damage the glass parts.
	Please connect the power supply according to the requirements on the name plate.
	Do not press the power cord. When there is a heavy thing on the power line or through the object, it may cause damage to the power cord and cause electric shock or fire.
	Please unplug the power plug from the socket to avoid long-term accumulation of dust if you do not use it for a long time, if the temperature is very high it may cause a fire.
	Do not pull the power cord strongly.
	Please disconnect the power supply and unplug the plug when moving it. Do not move it if there is power to avoid electric shock or fire.
	Please clean the glass parts before using it.
	When the glass device is connected, the interface should be uniformly applied to the vacuum grease to ensure the sealing effect.
	Please be careful when install the glass parts.
	Do not hit the glass.
	The viscosity of the circulating liquid must be less than 50cp when the reactor connected with dynamic temperature control system.
	Please start the circulating pump first, then start the Dynamic Temperature Control System.
Do not heating the reactor with steam.	
 <p>Warning</p>	Do not make the stirring paddle rotate at a high speed of no-load.
	Please contact us or searching for technical support when there's any abnormal situation.
	The maximum temperature difference between the jacket and the vessel is 80°C.
	Do not move the device when it is working.
	Be aware of scaling! Please wear glasses when dealing with the sample.
 <p>Warning</p>	Do not move it under high temperature!
	Do not refit this device.
	Do not use it outdoor. Get wet in the rain or splashing water on the surface may cause personnel injury.

4.2 Operating

Note

Note:

Before starting the stirring blade for the first time, pay attention to the relative position of the stirring paddle and the bottom of the vessel body to avoid collision.

- 1) Set the "power switch" on the rear panel of the control box to "1", the "speed display window" and the "temperature display window" of the control box are displaying the set stirring speed and the measured temperature in the vessel.

NOTE: Do not make the stirring paddle rotate at a high speed of no-load.

- 2) Add the solid material or liquid material to the vessel body slowly through corresponding feeding port.
- 3) Start rotation speed control through press 'Rotation' button on the control panel. And then adjust 'speed regulator' clockwise to make the stirring speed rise to the target value.
- 4) The reactor can be pumped to a negative pressure state according to some experimental requirements.
- 5) Start Dynamic Temperature Control System according as needed to achieve automatic temperature control.
- 6) After reaction completed, closed the Dynamic Temperature Control System when the material temperature in the vessel close to room temperature.

Note:

Do not close the control valve in the circulating pipe after close Dynamic Temperature Control System. Otherwise the vessel body will be broken when the temperature difference changes greatly.

Note

- 7) Power off the vacuum pump.
- 8) Turn the "speed regulator" counterclockwise to the left, press "Rotation" button, stop stirring, and set the power switch on the rear panel of the control box to "0".
- 9) Turn the black knob of the discharge valve (14) counterclockwise to discharge the material. After the discharge is completed, remove the discharge valve carefully and clean the reactor discharge valve and port. And then re-install the discharge valve.

Note:

- a) When discharging, rotate the black knob at the bottom of the discharge valve counterclockwise.
- b) After cleaning the glass, please use a clean soft cloth to wipe the liquid inside and outside the device to avoid contamination.

5. Maintenance and Management

Ensure to make the equipment running normally to increase its service life, users should do daily maintenance and management. Before maintenance, please disconnect equipment power supply, preparing tools, materials according to according to the specific requirements in the operation manual. Otherwise, it will cause electric shock or damage.

5.1 Management

Please follow following recommendations to keep the normal operation of the reactor:

- 1) Check the ambient temperature and humidity regularly to ensure a better working conditions.
- 2) Avoid severe vibration when moving it.
- 3) Beware of touch the vessel in case of broke glass parts.

If found the component aging or damage, please replace it timely according to the original specifications.

5.2 Cleaning

- 1) Before cleaning the unit, disconnect the power cord. Otherwise, it may cause electric shock or fire hazard.
- 2) Use soft cloth to clean the surface.
- 3) Cleaning of glass parts should be done according to laboratory regulations.
- 4) Beware of hazardous substance leaks on the surface of the vessel or enters it.
- 5) Do not use brushes, grinding powder, acid, petrol or other solvents to clean the surface of the body to avoid damage to the protective layer.
- 6) If the vessel and its connection glass parts' temperature is high, please disconnect the power until it is fully cooled.
- 7) In order to avoid damage of components, please do not press the monitor and knob forcedly when cleaning them.

5.3 Maintenance

The warranty of the equipment against manufacturing problems is of 12 months from the date of purchase. If the equipment has been used improperly, compensation for the damage caused will be charged.

Contact the Technical Service before obtaining professional maintenance to return the equipment to the factory:

1. Clean the machine so as not to damage the maintenance personnel.
2. Provide the description of the equipment failure.
3. Pack the machine properly to avoid damage during transport.


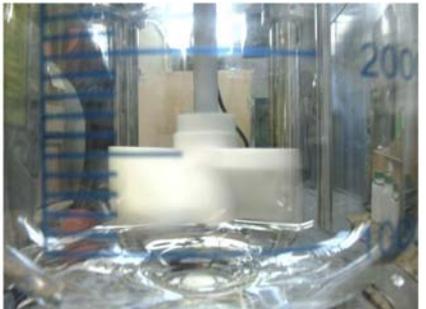
5.4 Troubleshooting

Table 5.3 Faults and troubleshooting

Faults	Reason	Troubleshooting	
No display on touch screen after power on	Power supply is not connected	Check the power supply circuit	
	* Fuse fault	Check the fuse and its installation	
	Power switch fault	Replace power switch	
	PCB board fault	Please contact us or professional maintenance personnel	
Motor rotate oppositely	Motor power line phase sequence fault	Change the line phase sequence	
Temperature display abnormal	PCB board fault	Please contact us or professional maintenance personnel	
	Improper wiring or poor contact		
	Sensor fault		
Rotation speed is '0', motor doesn't work	Motor fault		
	PCB board fault		
	Start button breakdown		
	Speed regulator breakdown		
Abnormal noise in stirring process	Stirring paddle installed improperly		Re-install stirring paddle
	Motor fault		Please contact us or professional maintenance personnel
	Debris in the vessel		Clean debris
Vacuum degree decrease	Sealing problem between glass parts	Re-install the glass parts	
	Sealing ring wear	Replace sealing ring	
	Sealing ring installed improperly	Re-install sealing ring	
	Vacuum hose aging	Replace vacuum hose	
	The vacuum capacity of the external vacuum pump is reduced	Replace vacuum pump	
Discharge valve leak air	When the material temperature in the vessel changes, the sealing of the discharge port and discharge valve will be affected due to different material thermal expansion coefficient difference	Regulate connecting flange	
	Sealing ring wear	Replace sealing ring	

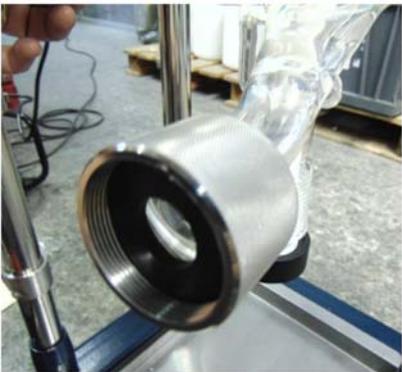Note: * The fuse is located below the power outlet of the rear panel of the controller.

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Coolant double effect DIN 29/32 L: 300mm

Addition funnel 250ml, compensated pressure DIN29 / 32

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Criothermostat “FRIGITERM-TFT-30”
Accessory for the coolant circuit
For regulable temperatures from -30 up to 100°C
Part number 3000547

Gable lid
Thermostatic bath accessory which limits evaporation
Part number 3001295

Thermostatic bath DIGITERM-TFT-200 -20
Accessory for the reactor
To regulate temperatures from ambient +5°C up to 200°C
Part number 3000550

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Water recirculator
"Interheater" Accessory of the
coolant circuit
To regulate temperature from +5° up to 90°C
Part number 6001491

Rotary vacuum pump
Flow volume 3.6 m³/h. Vacuum limit 0.1 mbar
Part number 5900620
spare oil for pump
Part number 5515030

Vacuum membrane pump N-820
Special for solvents
Flow volume 20 litres/minute. Vacuum limit 8 mbar. PTFE rotors. PTFE coated
membranes, valves in FFPM, motor protection level IP-44
Part number 5910009

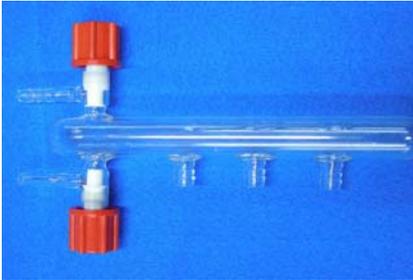
Scrubber unit
Glass trap in Pirex to avoid aspiration of liquids by means of the vacuum pump
Part number 4001611

Universal clamp
In dur-al with jaws in cork coated
Part number 7000403

Double nut
In chrome plated
Part number 7000408

Double glass valve
With three vacuum connection points (coolant, load & drip)
Part number 47078

Glass T
For vacuum connection
Part number 47077

Silicone pipe
External diameter 14 mm. Internal diameter 8 mm
Part number 46041

Corrugated tube
Reinforcement for the silicone tube
Part number 46079

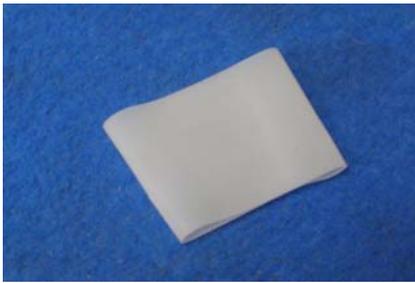
Fixing clamps
Bags of 100 units
Part number 28006 (unit)

PTFE bellows diameter 29/32 x 100mm
Part number 1001825

Conical connection cover 29/32 PTFE
Part number 28515

Glass sensor adapter for reactor
Combines with PH/T°C analyzer part number 4120800
& sensor part number 120801
Part number 4120802

Guide for PTFE stirrers 29/32 \varnothing 10 MM
High chemical resistance
Low vacuum (5mmHg)
Pressure (3-5 psi)
High performance PTFE/PEEK/glass precision bearing seal
Internal anti-whip sleeve
Part number 1001788

Addition funnel 250ml PTFE key Step 4mm
Part number 1001795

Load faucet 29/32 Step 4mm Olive 12/13
Part number 1001793

Stainless steel funnel with handle 200 mm
Part number 1000909

Support accessories trolley
Total height 90 cm
Lower platform dimensions: 70x70 cm
Upper platform dimensions: 70x40 cm
Part number 1001744 (unit)

Stainless steel bucket 15 litres
Part number 1001600

Large opening Hoffman clamp
Part number 1000064

50 litres reservoir
Part number 0106006

OTHER ACCESSORIES:

- 0046030 PTFE tube 10X8 (1 meter)
- 0046149 Silicon tube to unload 15x20. (1 meter)
- 0046150 Silvar tube 14.4x8 mm reinforced with fiberglass (5 meters)
- 0046156 Silvar tube 19x12 mm reinforced with fiberglass (5 meters)
- 0046084 Viton tube 10x14 mm (1 meter)
- 0046085 Viton tube 12x17 mm (1 meter)
- 0046041 Silicon tube 14x8 mm (5 meters)
- 9515031 Fluid 47 V50 4 litres (silicone oil)
- 1001853 PTFE corrugated tube DN12 connection Ø10 (1 meter)
- 1001915 High vacuum grease 50 gr.

7. Examples of Use of Support Trolleys

Frigiterm TFT & Interheater

Digiterm TFT & Interheater

Digiterm TFT, Vacuum pump, Scrubber & steel cubes
Vacuum pump, Scrubber and 50 litres reservoir

8. Examples of Application